
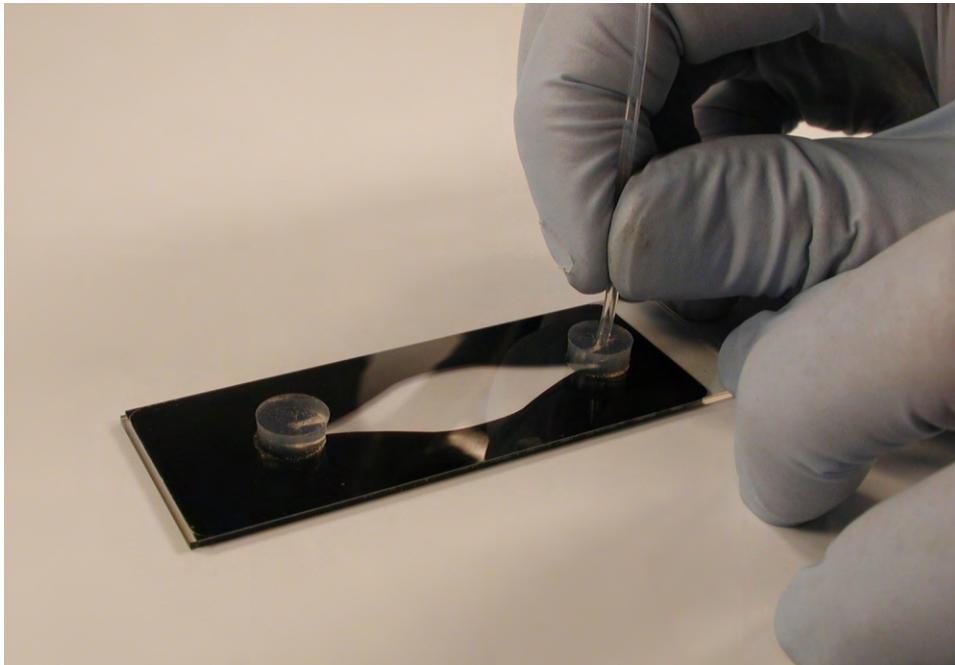
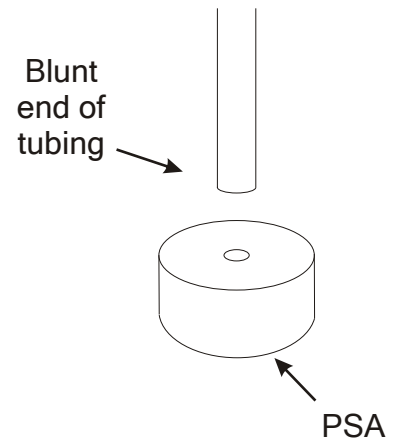


Press Fit Tubing Connectors
Item # 460003
20/PK

Incubation chambers and tubing
sold separately.

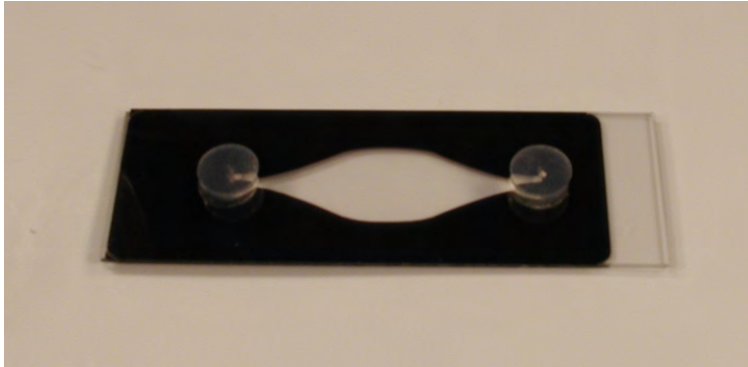

Remove paper liner from each connector.
Press one connector over each filling port
Trim tubing to desired length(s)
Press the tubing into the small hole in the top surface of each connector.
Tubing may be adapted to inlet connectors or syringes.
Recommended Tubing PTFE #24 (Cole Parmer, Inc. Item No. 06417-21)

Grace Bio-Labs, Inc.

Tel: 541-318-1208 / Fax: 541-318-0242
www.gracebio.com

