



IRIS9 Plus - Prestained Protein Ladder

24 DEC 2021

Catalog Number	Size	Range
PMI09-0500	500 µl	15 to 170 kDa

Storage Conditions

Stable for up to 2 weeks at 25°C.

Stable for up to 3 months at 4°C.

Stable for up to 24 months at -20°C.

Description

The IRIS9 Plus Prestained Protein Ladder is a combination of 9 pre-stained proteins with molecular weights from 15 to 170 kDa. The 9 recombinant proteins are covalently coupled with blue chromophore, while 1 green band at 15 kDa, a red band at 70 kDa and a newly designed peacock green band at 60 kDa serve as reference bands. The IRIS9 Plus Prestained Protein Ladder keeps track of the size and separation of proteins during SDS-polyacrylamide gel electrophoresis, approximating the target protein size and validating the Western transfer efficiency on PVDF, nylon, or nitrocellulose membranes.

Kit Content(s)

IRIS9 Plus Prestained Protein Ladder	500µl x 1 vial
--------------------------------------	----------------

Storage Buffer

0.2~0.4 mg/ml of each band of protein is stored in a mixture of solution that contains: 20 mM Trisphosphate at pH 7.5, 2% SDS, 0.2 mM Dithiothreitol, 3.6 M Urea, and 15% (v/v) Glycerol.

Required materials but not provided

- Vertical Electrophoresis system
- Power supplies
- Vortex or equivalent
- Microcentrifuge

Reaction Setup


3µl or 5µl per loading for clear visualization during electrophoresis on 15-well or 10-well mini-gel, respectively. 2.5µl per well for general Western transferring.





Guide for Molecular Weight Estimation (kDa)

Refer to the IRIS9 Plus Prestained Protein Ladder patterns in various electrophoresis conditions:

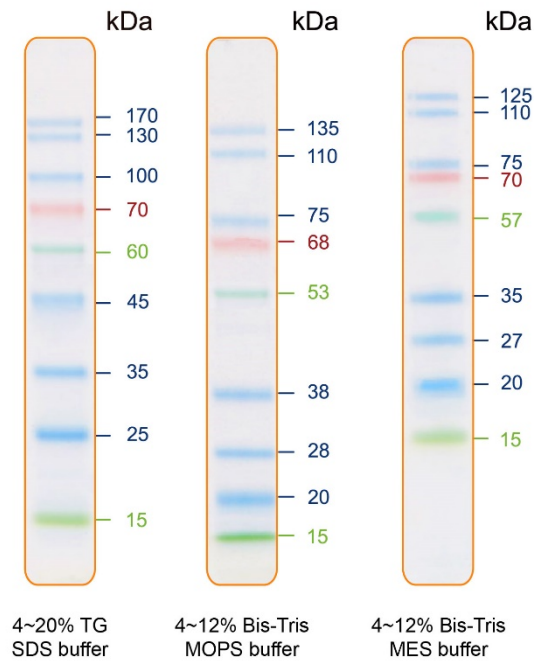

Important notes

The molecular weight of each protein (kDa) was measured against an unstained protein ladder in every electrophoresis condition. Additional data should be considered for a more accurate adjustment.