

# Carrez Clarification Reagent Kit

## Product Description

**Catalog #:** IDJ181-; 100 tests  
**Name :** **Carrez Clarification Reagent Kit**  
Kit content:  
Carrez Reagent I, 500 µl (Amber) #IJD18a  
Carrez Reagent II, 500 µl (Blue) #IJD18b

**Storage :** at Room Temperature (Z)



### I. Introduction:

Carrez clarification reagents are used to treat a wide variety of samples intended to be analyzed by enzymatic or other means. The reagents cause precipitation of proteins and fats, elimination of interference due to a number of redox compounds, which can affect assays and eliminates turbidity and emulsions particularly in food testing.

The treatment of samples with Carrez Clarification Reagents is quick, simple and dilutes samples by a minimal amount. The procedure is simpler, less expensive and takes less time than other methods such as spin filtering.

Most samples collected for analyses of small molecule analytes such as carbohydrates, alcohols, aldehydes and organic acids can be prepared using this reagent system.

Typical uses include glucose, fructose, sucrose, lactose, lactulose, maltose etc. ; ethanol, glycerol, sorbitol, acetaldehyde, lactate, nitrate.

However Carrez clarification is not suitable for samples in which enzymatic activities are to be quantified. These reagents must not be used to prepare samples for analytes such as ascorbate (vitamin C), citrate, urea (> ammonia), aconitate (> citrate).

### II. Application:

- Preparation of samples for further analysis (bioassay, HPLC, etc.)

### III. Sample Type:

- Foods (meats, milk and milk products, forage, cereals, etc.)
- Blood

### IV. Kit Contents

Carrez Reagent I, 500 µl (Amber) #IJD18a  
Carrez Reagent II, 500 µl (Blue) #IJD18b

### V. User supplied Reagents and Equipment

- Sample preparation equipment (homogenizer, Turrax, etc.)
- 0.22 µm syringe filters
- Buffer, pH 7.5-8.0 or 50 mM NaOH

FT-IDJ181

## VI. Storage and Handling:

Store kit at room temperature.

Avoid exposing Carrez Reagent I to strong acids as this can generate toxic cyanide gas.

Both reagents are supplied ready to use.

## VII. Protocol:

1. Take 10 mg of a sample and homogenize in 100  $\mu$ l of dH<sub>2</sub>O until a uniform suspension is formed. Transfer to a disposable microcentrifuge tube.

2. Add 5  $\mu$ l of Carrez Reagent I and vortex several sec. to mix.

Add 5  $\mu$ l Carrez Reagent II and vortex until homogeneous.

3. Centrifuge for 2 min. at 10,000 X g to pellet insolubles.

Transfer the clear supernatant to a fresh tube.

For analytes, which are present in relatively high abundance, dilute 100X with dH<sub>2</sub>O and use directly.

**Note:** For analytes present in small amounts where dilution of the sample may present detection problems, adjust the pH of the sample to pH 7.5-8.0 to precipitate excess Zinc ions from Reagent II. This can be accomplished with a buffer at a volume ratio of approximately 1:1 with sample. A brief centrifugation or filtration through a 0.22  $\mu$ m filter will remove the Zn(OH)<sub>2</sub> precipitate. Alternatively to minimize dilution, 50 mM NaOH can be used. Caution must be used in this case since the sample has a very low inherent buffer capacity and it is very easy to make the solution very alkaline, which can cause decomposition of many analytes.

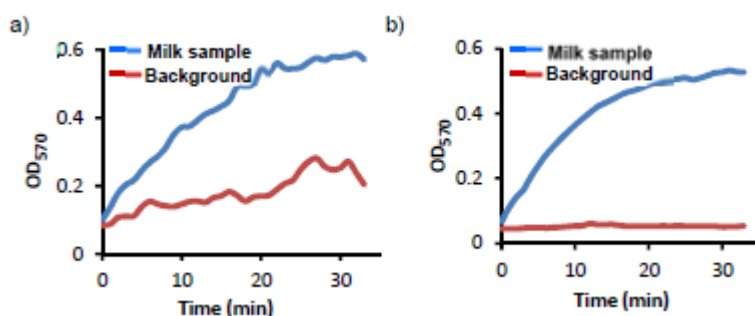


Figure: Comparison of analysis of glucose in milk using Carrez reagents. 100  $\mu$ l of fat-free milk untreated (a) or treated with 5  $\mu$ l of each Carrez Reagent sequentially (b) was centrifuged to remove precipitated material and diluted 100X. 10  $\mu$ l was added to a 96-well plate. Lactose in samples was hydrolyzed using  $\beta$ -galactosidase and glucose produced was analyzed using a Glucose Assay Kit ( ).

## VIII. RELATED PRODUCTS:

Deproteinizing Sample Preparation Kit ( ) 10K Spin Column ( )

## Ordering information

Catalog size quantities and prices may be found at <http://www.interchim.com>.

Please contact InterBioTech – Interchim for any other information

Hotline : +33(0)4 70 03 73 06 – [Interbiotech@interchim.com](mailto:Interbiotech@interchim.com)

**Disclaimer :** Materials are sold **for research use only**, and are not intended for food, drug, household, or cosmetic use. Interchim® is not liable for any damage resulting from handling or contact with this product.

Rev.Q04E