

# Product Information

## EvaEZ™ Fluorometric Polymerase Activity Assay Kit

**Catalog Number:** 29051

**Unit Size:** 2 x 1 mL (200 assays)

### Components

EvaEZ™ Fluorometric Polymerase Activity Mix contains primed substrate, EvaGreen® dye, dNTPs and ROX reference dye in polymerase buffer.

### Storage and Handling

Store at -20°C, protected from light. Product is stable for at least 6 months from date of receipt when stored as recommended.

### Product Description

EvaEZ™ Fluorometric Polymerase Activity Assay Kit provides an easy and accurate way to determine the activity of a nucleic acid polymerase without using radioisotopes. It contains EvaGreen® dye together with a primed template, dNTPs, MgCl<sub>2</sub>, and ROX reference dye (high ROX) in a Tris buffer system. In the presence of DNA polymerase activity, the primer will be extended to form a double stranded product that can bind EvaGreen dye, resulting in an increase of fluorescence (Figure 1). The rate of increase of fluorescence is positively correlated to the activity of polymerase (see example in Protocol B). The assay was developed for the measurement of Taq DNA polymerase activity. It can also be used for other DNA polymerases, such as Pfu, Vent, Phusion®, Bst, Phi29, MMLV, AMV, SuperScript®, T4 DNA polymerase, T7 DNA polymerase, Klenow, and E. coli DNA polymerase I. The activity assay can be conducted at temperatures from 4°C to 75°C.

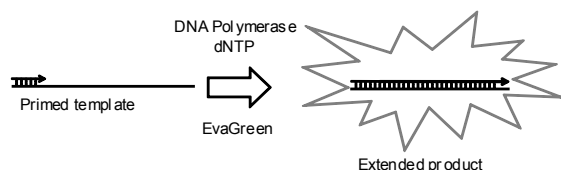


Figure 1. Schematic overview of EvaEZ fluorometric polymerase activity assay.

### Recommended Protocols

For researchers with basic knowledge of enzyme kinetics and activity assays. Protocol A describes how to measure fluorescence change by polymerization. Protocols B and C describe two different methods for using the fluorescence change by polymerization to calculate units of polymerase activity.

**Protocol A: Measurement of Fluorescence Change by Polymerization.**

- Combine the following components in a real-time qPCR reaction tube on ice:  
10 µl 2X EvaEZ Polymerase Activity Mix  
9 µl H<sub>2</sub>O  
1 µl DNA polymerase sample  
Gently mix the reaction components well.

- Quickly place the reaction tube into a real time qPCR instrument. Run an isothermal program at the designated temperature for the polymerase being tested (e.g. 37 °C for Klenow) for 60 minutes. Measure fluorescence in Channel 1 (i.e. the channel for FAM or EvaGreen).
- Line A in Figure 2 is a typical reaction trace recorded by a real time qPCR instrument (e.g. ABI 7900). The X-axis is time in minutes; the Y-axis is fluorescence. Make a contingent (Line B) that passes through time 0. The initial rate of fluorescence change (fluorescence unit/minute) resulting from the polymerase activity is represented by the slope of Line B.

**Note 1:** Depending on the instrument used, there may be a lag between the starting time of polymerization and the starting time when the fluorescence is collected.

**Note 2:** Reactions can also be carried out in a fluorometer cuvette or a plate reader where the temperature can be accurately controlled.

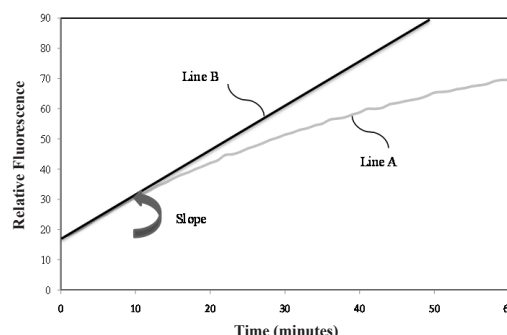


Figure 2. A typical reaction profile of EvaEZ fluorometric polymerase activity assay.

**Protocol B: Activity Determination of an Unknown Polymerase Sample Based on a Polymerase Standard.**

- Make a serial dilution of a standard DNA polymerase of known activity.
- Monitor the fluorescence of each dilution using Protocol A. Figure 3 shows a typical collection of reaction curves from a dilution series.
- Determine the initial rate of fluorescence increase for each polymerase dilution by calculating the slope of the initial linear portion of the curve before fluorescence begins to plateau (see Protocol A, step 3 and Figure 2).
- Plot the initial slope for each standard dilution against the unit activity of the standard DNA polymerase to generate a standard curve. A Taq polymerase standard curve is shown in Figure 4.
- Use the linear range of the plot from step 4 as a standard curve. Measure the initial rate of an unknown polymerase sample using Protocol A.
  - If the rate falls in the linear range of the standard curve, use the standard curve to calculate the activity of the unknown polymerase sample.
  - If the rate is outside the linear range, adjust the concentration of the unknown polymerase sample so that the initial rate falls into the linear range of the standard curve. Factor in the dilution factor when calculating the activity.

Continued next page

Protocol C: Activity Determination of an Unknown Polymerase Sample Based on the Number of Nucleotides Synthesized.

1. Make a serial dilution of an unknown DNA polymerase and measure the initial rates according to Protocol A.
2. Select an enzyme concentration from the dilution series that is in the linear response region and determine the slope of the initial fluorescence change as described in Protocol A. By definition, as shown in Figure 5:

$$\Delta F = \text{slope (fluorescence unit/min)} * 60 \text{ (min)}$$

3. Separately, run polymerization reactions with enzyme of a saturating concentration and a no enzyme control respectively according to the procedure described in Protocol A.
4. Derive the maximum fluorescence change ( $\Delta F_{\text{max}}$ ) by subtracting the fluorescence at the 60th minute of the no enzyme control from the fluorescence at the 60th minute of the saturating enzyme.

The number of nucleotides synthesized by the polymerase at the selected concentration in 60 minutes under kit conditions is:

$$(\Delta F / \Delta F_{\text{max}}) * 270 \text{ pmole}$$

Note: 270 pmol is the the number of nucleotides synthesized in a 20  $\mu\text{L}$  EvaEZ reaction with saturated enzyme.

References:

1. Irwin H. Segel, I. H. Segel. Enzyme Kinetics: Behavior and Analysis of Rapid Equilibrium and Steady-State Enzyme Systems. 1976.
2. Mao F, Leung WY, Xin X. (2007). Characterization of EvaGreen and the implication of its physicochemical properties for qPCR applications. BMC Biotechnol. 7, 76.

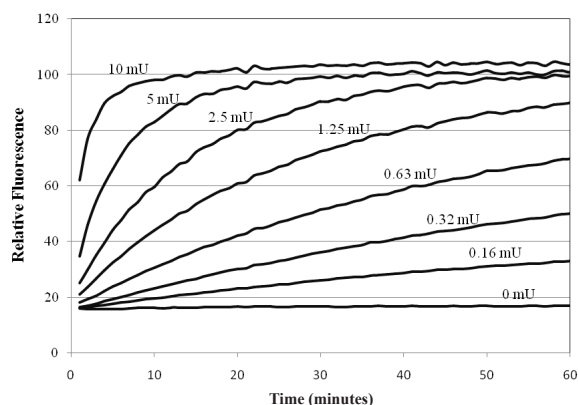


Figure 3. Typical reaction curves from a polymerase dilution series.

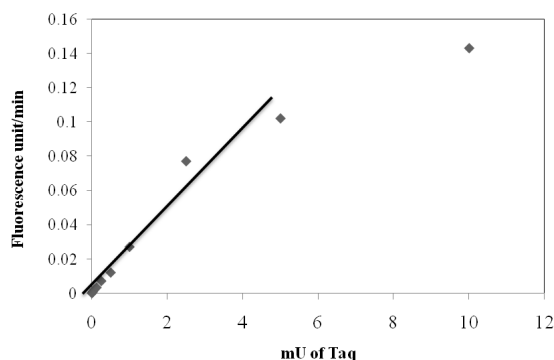


Figure 4. Titration of Taq DNA Polymerase Activity.

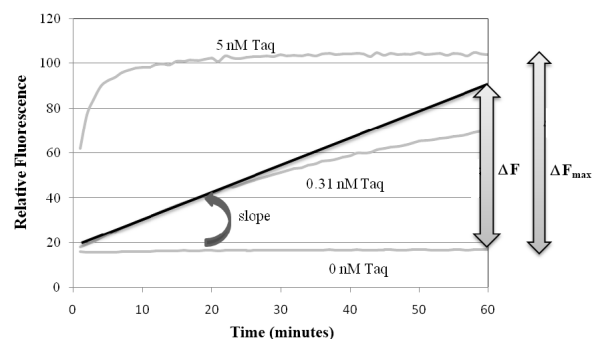


Figure 5. Activity determination based on the number of nucleotides synthesized.

Related Products

Catalog number	Product	
31000-T	EvaGreen dye, 20X in water, trial size, 1 mL	1 mL
29050	Cheetah HotStart Taq DNA polymerase, 500 U	500U
31020-T	Fast Plus EvaGreen Master Mix, no ROX, 100 rxn	2x1ml
31014-T	Fast Plus EvaGreen Master Mix, low ROX, 100 rxn	2x1ml
31015-T	Fast Plus EvaGreen Master Mix, high ROX, 100 rxn	2x1ml
29054	HotStart Polymerase Modification Kit	1 kit
29053	WarmStart Polymerase Modification Kit	1 kit
41003	GelRed Nucleic Acid Gel Stain, 10,000X in H <sub>2</sub> O	0.5 mL
41005	GelGreen Nucleic Acid Gel Stain, 10,000X in H <sub>2</sub> O	0.5 mL

Please visit our website at [www.biotium.com](http://www.biotium.com) for information on our life science research products, including environmentally friendly EvaGreen® qPCR master mixes, fluorescent CF™ dye antibody conjugates and reactive dyes, apoptosis reagents, fluorescent probes, and kits for cell biology research.

Materials from Biotium are sold for research use only, and are not intended for food, drug, household, or cosmetic use.

EvaEZ™ is covered under pending US and international patents. EvaGreen® dye and applications are covered under US patent nos. 7,803,943 and 7,776,567 and pending international patents. Our products are not available for resale or other commercial uses without a specific agreement from Biotium, Inc. We welcome inquiries about licensing the use of our dyes, trademarks or technologies. Please submit inquiries by e-mail to [btinfo@biotium.com](mailto:btinfo@biotium.com). EvaGreen® is a registered trademark of Biotium, Inc. Phusion® is a registered trademark of Thermo Fisher Scientific. SuperScript® is a registered trademark of Invitrogen.