



ISO 9001:2000 CERTIFIED
ISO 13485:2003 CERTIFIED



Freezing Media, Bacterial

N400-25PK

Directions for Use

Components: 0.5 ml aliquot / cryotube

- LB broth (Miller)
- 30 % glycerol

Procedure:

1. Add 1 ml of saturated culture to the vial.
2. Vortex well.
3. Store -70 °C.
4. To inoculate fresh culture, soften briefly on ice, remove small chunk and add to fresh media.

Storage/Stability:

Store at 2-8 °C.



211 Avenue J.F. Kennedy - BP 1140

03103 Montluçon cedex - France

Tél 33 (0)4 70 03 88 55 - Fax 33 (0)4 70 03 82 60

e-mail interchim@interchim.com - www.interchim.com