



Quality & Excellence since 1996

FD Rapid GolgiStain™ Kit

A complete Golgi-Cox staining system for the study of the morphology of neurons and glia

User Manual

PK 401/401A, Version 2012-02

FOR *IN VITRO* RESEARCH USE ONLY
not for diagnostic or other uses

FD NeuroTechnologies
Consulting & Services, Inc.

© 2002-2012 All Rights Reserved

Contents

I.	Introduction	4
II.	Kit Contents	5
III.	Materials Required but Not Included	5
IV.	Safety and Handling Precautions	6
V.	Tissue Preparation	6
VI.	Staining Procedure	11
VII.	References	13

I. Introduction

Golgi-Cox impregnation^{2,5} has been one of the most effective techniques for studying both the normal and abnormal morphology of neurons as well as glia. Using the Golgi technique, subtle morphological alterations in neuronal dendrites and dendritic spines have been discovered in the brains of animals treated with drugs as well as in the postmortem brains of patients with neurological diseases^{1,3}. However, the reliability and time-consuming process of Golgi staining have been major obstacles to the widespread application of this technique.

FD Rapid GolgiStain™ kit is designed based on the principle of the methods described by Ramón-Moliner⁵ and Glaser and Van der Loos⁴. This kit has not only dramatically improved and simplified the Golgi-Cox technique but has also proven to be extremely reliable and sensitive for demonstrating morphological details of neurons and glia, especially dendritic spines. The FD Rapid GolgiStain™ kit has been tested extensively on the brains from several species of animals as well as on the specimens of postmortem human brains (for photo samples and references using this kit, please visit our web site at www.fdtype.com).

II. Kit Contents

Store at room temperature

	<u>PK401A</u>	<u>PK401</u>
Solution A	125 ml	250 ml
Solution B	125 ml	250 ml
Solution C	125 ml x 2	250 ml x 2
Solution D	125 ml	250 ml
Solution E	125 ml	250 ml
Glass specimen retriever	2	2
Natural hair paintbrush	2	2
Dropping bottle	1	1
Plastic forceps	1	1
User manual	1	1

III. Materials Required but Not Included

1. Double distilled or Milli-Q water
2. Plastic/glass tubes or vials
3. Histological supplies and equipment:
 - Gelatin-coated microscope slides
 - Coverslips
 - Staining jars
 - Ethanol
 - Xylene
 - Permount[®]
 - A light microscope.

Permount[®] is a registered trademark of Fisher Scientific.

IV. Safety and Handling Precautions*

1. FD Rapid GolgiStain™ kit is made for *in vitro* research use only and not for drug, diagnostic or other uses.
2. The kit contains reagents that are toxic and harmful in contact with skin or by inhalation and may be fatal if ingested. Do not pipette by mouth. Avoid inhalation and contact with skin and eyes. In case of contact, wash immediately with generous amounts of water and seek medical advice. If swallowed, wash out mouth with water and immediately call a physician.
3. Perform experiment under a chemical hood. **Wear suitable protective clothing, gloves and eye/face protection while handling kit reagents.** Wash hands thoroughly after performing the experiment.

*Material safety data sheet (MSDS) is available at www.fdneurotech.com.

V. Tissue Preparation

The following instructions must be read before using this kit.

- All containers (plastic preferred) to be used should be cleansed and rinsed with distilled water.
- Do not use metal implements whenever Solutions A and B are present.
- Keep containers tightly closed at all times.
- Tissues treated with Solutions A and B, including sections should be protected from light whenever possible.

- The following procedure should be performed at room temperature unless specifically indicated.

 Note

FD Rapid GolgiStain™ kit has been proven to produce the best results in animal and postmortem human brain tissues prepared according to the following procedures. However, it may be used in tissue specimens prepared differently (please contact FD NeuroTechnologies for more information before use).

1. Experimental animals should be deeply anesthetized before killing. The animal brain (or postmortem human specimens) should be removed from the skull as quickly as possible but handled carefully to avoid damage or pressing of the tissue.

 Note

- *Perfusion of animals is not necessary. However, if you do perfuse animals (e.g. with 4% paraformaldehyde for less than 5 minutes), tissue must not be postfixed. Important!*
 - *Large brain specimens should be sliced with a sharp blade into blocks of approximately 10 mm thickness.*
2. Rinse tissue briefly in double distilled or Milli-Q water to remove blood from the surface.
 3. Immerse tissue in the impregnation solution, made by mixing equal volumes of Solutions A and B, and store at room temperature for 2 weeks in the dark*. Replace the impregnation solution after the first 6 hours of immersion or on the next day.

** A 2-week impregnation time is satisfactory in most cases. However, variations in type and actual size of tissue may require a shorter or longer duration of impregnation to obtain the best results. The ideal time should be obtained by trial for each type of tissue, but 3 weeks is often sufficient for most tissues. Note that prolonging the impregnation time may increase background staining.*

Note

- *For the best results, the mixture of Solution A and B should be prepared at least 24 hours prior to use and left unstirred.*
- *It is important to use the top part of solution that is free of precipitate.*
- *The impregnation solution may be stored at room temperature for up to 1 month in the dark before use.*
- *Use at least 5 ml of the impregnation solution for each cubic cm of the tissue to be studied (i.e., the volume of the impregnation solution should be at least five times that of the tissue). Note that use a lesser volume of impregnation solution may decrease the sensitivity and reliability of staining.*

Warning:

Solutions A and B (containing mercuric chloride, potassium dichromate and potassium chromate) are toxic in contact with skin and may be fatal if swallowed. The experiment should be performed under a chemical hood. Wear suitable protective clothing, gloves and eye/face protection while handling the reagents. DO NOT POUR THE WASTE OF SOLUTIONS A AND B INTO THE

SINK. Collect the waste of these solutions in a bottle and call your safety office or a licensed professional waste disposal service to dispose of this material.

4. Transfer tissue into Solution C and store at 4°C in the dark for at least 48 hours (up to 1 week). Replace the solution at least once after the first 24 hours of immersion or on the next day.
5. Sections (80 to 240 μm in thickness) can be best cut on a cryostat at -20°C to -22°C. Other types of microtomes, including sliding microtome and vibratome, can also be used (for tips, cf. pages 10 & 11). Each section should then be transferred with a glass specimen retriever (provided) and mounted on gelatin-coated microscope slides with Solution C (a dropping bottle is provided for easy dropping of Solution C onto the slides). Excess solution left on slide should be suctioned with a Pasteur pipette and then absorbed with a strip of filter paper (solution on slide must be wiped away as much as possible or sections will fall off slides). Allow sections to dry naturally at room temperature (do not use a fan or hot plate). For the best results, sections should be processed as soon as possible, but may be stored in a slide box at room temperature in the dark for up to 3 days.

 Note

- *To prevent tissue from ice crystal damage and to preserve the best possible cell morphology, tissue should*

be frozen rapidly before sectioning with a cryostat. For example, tissue may be rapidly frozen as described below: place tissue in a plastic spoon and slowly dip into iso-pentane pre-cooled with dry ice (for the best results, temperature should be kept below -70°C and the dipping should take about 1 min, the slower the better). After the tissue is completely immersed in iso-pentane, keep it in iso-pentane for a few seconds and then place it on dry ice for another minute to ensure that the tissue is well frozen. Do not let tissue thaw before sections are cut.

- The types of cryostat may vary, but all types should be able to cut thick sections (e.g. $100\ \mu\text{m}$). Contact FD Neuro-Technologies for technical assistance.
- If the cryostat has only one temperature control, set the cryostat temperature to -22°C at least 4 hours before cutting. If the cryostat has 2 temperature settings, set the chamber temperature 1 degree lower (colder) than that of the specimen head. Please note that -22°C is satisfactory in most cases. However, variations in type of cryostat and tissue may require a higher or lower chamber temperature in order to cut sections smoothly and without shattering.
- The following tips should be helpful for sectioning with a cryostat: 1) Mount tissue on specimen disc (also called chuck or stage) with distilled water, but make sure that the tissue is not thawed (may be done on dry ice). Tissue may also be mounted with any type of tissue freezing medium, including OCT, but avoid cutting through the medium. Do not embed the tissue in OCT. If the tissue has to be embedded for cutting, use TFM (TBS, Durham, NC, USA, Cat. #TFM-5) instead; 2) After the tissue is mounted on specimen disc, place the tissue on dry ice for 10 minutes, and then immediately install the specimen disc together with

the tissue on a cryostat and wait 5 minutes before cutting;
3) Cut a few sections (do not use the anti-roll plate). If the tissue is too cold, e.g. sections show cracks which are parallel to the blade, wait a few more minutes before the next trial. Otherwise continue to cut.

- *The impregnated brain may also be embedded in agarose or gelatin and cut with a vibratome. However, for collecting sections (filling the chamber), Solution C must be used. Otherwise, sections may crack upon drying.*
- *To cut the impregnated tissue with a sliding microtome, both the stage and knife (or blade) need to be maintained at low temperature (below 0°C). It is also important that sections should be mounted with solution C as described in the User Manual.*

VI. Staining Procedure

Do not let sections dry out between any steps during the staining and before coverslipping.

1. Rinse sections in double distilled or Milli-Q water 2 times, 2 minutes each.
2. Place sections in a mixture consisting of 1 part **Solution D**, 1 part **Solution E** and 2 parts double distilled or Milli-Q water for 10 minutes.

e.g. Solution D	10 ml
Solution E	10 ml
Double distilled water	20 ml

 Note

- *The working solution should be prepared just before use and may be used for up to 200 sections (e.g. mouse brain) per 100 ml, depending on the size of sections.*
 - *The bottle and staining jar containing the working solution must be covered to prevent vaporization of the reagent.*
 - *For the best results, the working solution should be stirred frequently during incubation.*
3. Rinse sections in double distilled or Milli-Q water 2 times, 4 minutes each (distilled water should be renewed frequently).
 4. Counterstain sections with cresyl violet or thionin (optional step).

 Note

- *For counterstaining, step 3 must be prolonged, e.g. instead of 2 times and 4 minutes each, wash 4 times and 5 minutes each or longer.*
5. Dehydrate sections in 50%, 75% and 95% ethanol, 4 minutes each (do not skip any step).
 6. Dehydrate sections in absolute ethanol, 4 times, 4 minutes each (do not prolong).
 7. Clear in xylene, 3 times, 4 minutes each, and coverslip with Permount[®].

 Note

- *Sections may be temporally stored in xylene for a few*

hours before coverslipping.

- *For the best results, use undiluted PermOUNT[®].*
- *Golgi-stained sections should be protected from light whenever possible.*

VII. References

1. Robinson TE and Kolb B. Persistent structural modification in nucleus accumbens and prefrontal cortex neurons produced by previous experience with amphetamine. *J. Neurosci.* 17:8491-8497, 1997.
2. Corsi P. Camillo Golgi's morphological approach to neuroanatomy. In Masland RL, Portera-Sanchez A and Toffano G (eds.), *Neuroplasticity: a new therapeutic tool in the CNS pathology*, pp 1-7. Berlin: Springer, 1987.
3. Graveland GA, Williams RS and DiFiglia M. Evidence for degenerative and regenerative changes in neostriatal spiny neurons in Huntington's disease. *Science* 227:770-773, 1985.
4. Glaser ME and Van der Loos H. Analysis of thick brain sections by obverse-reverse computer microscopy: application of a new, high clarity Golgi-Nissl stain. *J. Neurosci. Methods* 4:117-125, 1981.
5. Ramón-Moliner E. The Golgi-Cox technique. In Nauta WJH and Ebbesson SOE (eds.), *Contemporary Methods in Neuroanatomy*. pp 32-55, New York: Springer, 1970.

References using this kit:

1. Adlard PA, Bica L, White AR, Nurjono M, Filiz G, Crouch PJ, Donnelly PS, Cappai R, Finkelstein DI, Bush AI. Metal ionophore treatment restores dendritic spine density and synaptic protein levels in a mouse model of Alzheimer's disease. **PLoS One.** 6(3):e17669, 2011.
2. Alberi L, Liu S, Wang Y, Badie R, Smith-Hicks C, Wu J, Pierfelice TJ, Abazyan B, Mattson MP, Kuhl D, Pletnikov M, Worley PF, Gaiano N. Activity-induced Notch signaling in neurons requires Arc/Arg3.1 and is essential for synaptic plasticity in hippocampal networks. **Neuron.** 69(3):437-44, 2011.
3. Anderson-White BR, Ivey FD, Cheng K, Szatanek T, Lorestani A, Beckers CJ, Ferguson DJP, Sahoo N, and Gubbels M-J. A family of intermediate filament-like proteins is sequentially assembled into the cytoskeleton of *Toxoplasma gondii*. **Cell. Microbiol.** 13(1):18-31, 2011.
4. Andres RH, Horie N, Slikker W, Keren-Gill H, Zhan K, Sun G, Manley NC, Pereira MP, Sheikh LA, McMillan EL, Schaar BT, Svendsen CN,

- Bliss TM, and Steinberg GK. Human neural stem cells enhance structural plasticity and axonal transport in the ischaemic brain. **Brain**. 134(6):1777-89, 2011.
5. Antzoulatos E, Jakowec MW, Petzinger GM, and Wood RI. MPTP neurotoxicity and testosterone induce dendritic remodeling of striatal medium spiny neurons in the C57Bl/6 mouse. **Parkinson's Dis**. doi:10.4061/2011/138471, 2011.
 6. Antzoulatos E, Jakowec MW, Petzinger GM, Wood RI. MPTP Neurotoxicity and Testosterone Induce Dendritic Remodeling of Striatal Medium Spiny Neurons in the C57Bl/6 Mouse. **Parkinsons Dis**. 2011:138471, 2011.
 7. Ayhan Y, Abazyan B, Nomura J, Kim R, Ladenheim B, Krasnova IN, Sawa A, Margolis RL, Cadet JL, Mori S, Vogel MW, Ross CA, and Pletnikov MV. Differential effects of prenatal and postnatal expressions of mutant human DISC1 on neurobehavioral phenotypes in transgenic mice: evidence for neurodevelopmental origin of major psychiatric disorders. **Mol. Psychiat**. 16(3):293-306, 2011.
 8. Babus LW, Little EM, Keenoy KE, Minami SS, Chen E, Song JM, Caviness J, Koo SY, Pak DT, Rebeck GW, Turner RS, Hoe HS. Decreased dendritic spine density and abnormal spine morphology in Fyn knockout mice. **Brain Res**. 1415:96-102, 2011.
 9. Bedrosian TA, Fonken LK, Walton JC, Haim A, Nelson RJ. Dim light at night provokes depression-like behaviors and reduces CA1 dendritic spine density in female hamsters. **Psychoneuroendocrinology**. 36(7):1062-9, 2011.
 10. Blugeot A, Rivat C, Bouvier E, Molet J, Mouchard A, Zeau B, Bernard C, Benoliel JJ, Becker C. Vulnerability to depression: from brain neuroplasticity to identification of biomarkers. **J. Neurosci**. 31(36):12889-99, 2011.
 11. Bongmba OYN, Martinez LA, Elhardt ME, Butler K, and Tejada-Simon MV. Modulation of dendritic spines and synaptic function by Rac1: a possible link to Fragile X syndrome pathology. **Brain Res**. 1399:79-95, 2011.
 12. Borgesius NZ, van Woerden GM, Buitendijk GH, Keijzer N, Jaarsma D, Hoogenraad CC, Elgersma Y. β CaMKII plays a nonenzymatic role in hippocampal synaptic plasticity and learning by targeting α CaMKII to synapses. **J. Neurosci**. 31(28):10141-8, 2011.
 13. Büdefeld T, Tobet SA, Majdic G. Altered position of cell bodies and fibers in the ventromedial region in SF-1 knockout mice. **Exp Neurol**. 232(2):176-84, 2011.
 14. Caito SW, Milatovic D, Hill KE, Aschner M, Burk RF, and Valentine WM. Progression of neurodegeneration and morphologic changes in the brains of juvenile mice with selenoprotein P deleted. **Brain Res**. 1398:1-12, 2011.
 15. Chen CC, Tam D, Brumberg JC. Sensory deprivation differentially impacts the dendritic development of pyramidal versus non-pyramidal neurons in layer 6 of mouse barrel cortex. **Brain Struct Funct**. [Epub ahead

of print] PMID: 21861159, 2011.

16. Chen C-Y, Lin C-W, Chang C-Y, Jiang S-T, and Hsueh Y-P. Sarm1, a negative regulator of innate immunity, interacts with syndecan-2 and regulates neuronal morphology. **J. Cell Biol.** 193(4):769-84, 2011.
17. Crupi R, Mazzon E, Marino A, La Spada G, Bramanti P, Battaglia F, Cuzzocrea S, and Spina E. Hypericum perforatum treatment: effect on behavior and neurogenesis in a chronic stress model in mice. **BMC Comp. Alt. Med.** 11:7, 2011.
18. DeVito LM, Balu DT, Kanter BR, Lykken C, Basu AC, Coyle JT, and Eichenbaum H. Serine racemase deletion disrupts memory for order and alters cortical dendritic morphology. **Genes, Brain, & Behav.** 10(2):210-22, 2011.
19. Dumanis SB, Cha H-J, Song JM, Trotter JH, Spitzer M, Lee J-Y, Weeber EJ, Turner RS, Pak DTS, Rebeck GW, and Hoe H-S. ApoE receptor 2 regulates synapse and dendritic spine formation. **PLoS One.** 6(2):e17203, 2011.
20. Duque A, Coman D, Carlyle BC, Bordner KA, George ED, Papademetris X, Hyder F, Simen AA. Neuroanatomical changes in a mouse model of early life neglect. **Brain Struct Funct.** [Epub ahead of print] PMID: 21984312, 2011.
21. Eiland L, Ramroop J, Hill MN, Manley J, McEwen BS. Chronic juvenile stress produces corticolimbic dendritic architectural remodeling and modulates emotional behavior in male and female rats. **Psychoneuroendocrinology.** [Epub ahead of print] PMID: 21658845, 2011.
22. Gao C, Frausto SF, Guedea AL, Tronson NC, Jovasevic V, Leaderbrand K, Corcoran KA, Guzmán YF, Swanson GT, and Radulovic J. IQGAP1 regulates NR2A signaling, spine density, and cognitive processes. **J. Neurosci.** 31(23):8533-42, 2011.
23. Gao X, and Chen J. Mild traumatic brain injury results in extensive neuronal degeneration in the cerebral cortex. **J. Neuropathol. & Exp. Neurol.** 70(3):183-91, 2011.
24. Gao X, Deng P, Xu ZC, Chen J. Moderate traumatic brain injury causes acute dendritic and synaptic degeneration in the hippocampal dentate gyrus. **PLoS One.** 6(9):e24566, 2011.
25. Hill MN, Hillard CJ, and McEwen BS. Alterations in corticolimbic dendritic morphology and emotional behavior in cannabinoid CB₁ receptor-deficient mice parallel the effects of chronic stress. **Cereb. Cortex.** 21(9):2056-64, 2011.
26. Hoffman AN, Krigbaum A, Ortiz JB, Mika A, Hutchinson KM, Bimonte-Nelson HA, Conrad CD. Recovery after chronic stress within spatial reference and working memory domains: correspondence with hippocampal morphology. **Eur J Neurosci.** 34(6):1023-30, 2011.
27. Huang W, Zhou Z, Asrar S, Henkelman M, Xie W, and Jia Z. p21-activated kinases 1 and 3 control brain size through coordinating neuronal complexity and synaptic properties. **Mol. and Cell Biol.** 31(3):388-403, 2011.

28. Jespersen S, Leigland L, Cornea A, Kroenke C. Determination of Axonal and Dendritic Orientation Distributions within the Developing Cerebral Cortex by Diffusion Tensor Imaging. **IEEE Trans Med Imaging**. [Epub ahead of print] PMID: 21768045, 2011.
29. Kambe T, Motoi Y, Inoue R, Kojima N, Tada N, Kimura T, Sahara N, Yamashita S, Mizoroki T, Takashima A, Shimada K, Ishiguro K, Mizuma H, Onoe H, Mizuno Y, and Hattori N. Differential regional distribution of phosphorylated tau and synapse loss in the nucleus accumbens in tauopathy model mice. **Neurobiol. Dis.** 42(3):404-14, 2011.
30. Kucukdereli H, Allen NJ, Lee AT, Feng A, Ozlu MI, Conatser LM, Chakraborty C, Workman G, Weaver M, Sage EH, Barres BA, Eroglu C. Control of excitatory CNS synaptogenesis by astrocyte-secreted proteins Hevin and SPARC. **Proc Natl Acad Sci U S A.** 108(32):E440-9, 2011.
31. Lancaster MA, Gopal DJ, Kim J, Saleem SN, Silhavy JL, Louie CM, Thacker BE, Williams Y, Zaki MS, Gleeson JG. Defective Wnt-dependent cerebellar midline fusion in a mouse model of Joubert syndrome. **Nature Med.** 17(6):726-31, 2011.
32. Lawrence RC, Otero NK, Kelly SJ. Selective effects of perinatal ethanol exposure in medial prefrontal cortex and nucleus accumbens. **Neurotoxicol Teratol.** [Epub ahead of print] PMID: 21871563, 2011.
33. Lioy DT, Garg SK, Monaghan CE, Raber J, Foust KD, Kaspar BK, Hirrlinger PG, Kirchhoff F, Bissonnette JM, Ballas N, Mandel G. A role for glia in the progression of Rett's syndrome. **Nature.** 475(7357):497-500, 2011.
34. Luine VN, Wallace ME, Frankfurt M. Age-related deficits in spatial memory and hippocampal spines in virgin, female Fischer 344 rats. **Curr Gerontol Geriatr Res.** 2011:316386, 2011.
35. Martin BJ, Naughton BJ, Thirtamara-Rajamani K, Yoon DJ, Han DD, Devries AC, Gu HH. Dopamine transporter inhibition is necessary for cocaine-induced increases in dendritic spine density in the nucleus accumbens. **Synapse.** 65(6):490-6, 2011.
36. Matrisciano F, Busceti CL, Bucci D, Orlando R, Caruso A, Molinaro G, Cappuccio I, Rizzo B, Gradini R, Motolese M, Caraci F, Copani A, Scaccianoce S, Melchiorri D, Bruno V, Battaglia G, and Nicoletti F. Induction of the Wnt antagonist Dickkopf-1 is involved in stress-induced hippocampal damage. **PLoS One.** 6(1):e16447, 2011.
37. Odajima J, Wills ZP, Ndassa YM, Terunuma M, Kretschmannova K, Deeb TZ, Geng Y, Gawrzak S, Quadros IM, Newman J, Das M, Jecrois ME, Yu Q, Li N, Bienvenu F, Moss SJ, Greenberg ME, Marto JA, Sicinski P. Cyclin E Constrains Cdk5 Activity to Regulate Synaptic Plasticity and Memory Formation. **Dev Cell.** [Epub ahead of print] PMID: 21944720, 2011.
38. Onksen JL, Brown EJ, and Blendy JA. Selective deletion of a cell cycle checkpoint kinase (ATR) reduces neurogenesis and alters responses in rodent models of behavioral affect. **Neuropsychopharmacol.** 36:960-9, 2011.

39. Paterniti I, Mazzon E, Gil C, Impellizzeri D, Palomo V, Redondo M, Perez DI, Esposito E, Martinez A, and Cuzzocrea S. PDE7 inhibitors: new potential drugs for the therapy of spinal cord injury. **PLoS One**. 6(1):e15937, 2011.
40. Perez-cruz C, Nolte MW, van Gaalen MM, Rustay NR, Termont A, Tanghe A, Kirchhoff F, and Ebert U. Reduced spine density in specific regions of CA1 pyramidal neurons in two transgenic mouse models of Alzheimer's disease. **J. Neurosci**. 31(10):3926-34, 2011.
41. Qin M, Entezam A, Usdin K, Huang T, Liu Z-H, Hoffman GE, and Smith CB. A mouse model of the fragile X permutation: effects on behavior, dendrite morphology, and regional rates of cerebral protein synthesis. **Neurobiol. Dis**. 42(1):85-98, 2011.
42. Qin M, Xia Z, Huang T, Smith CB. Effects of chronic immobilization stress on anxiety-like behavior and basolateral amygdala morphology in Fmr1 knockout mice. **Neuroscience**. 194:282-90, 2011.
43. Raman D, Milatovic SZ, Milatovic D, Splittgerber R, Fan GH, Richmond A. Chemokines, macrophage inflammatory protein-2 and stromal cell-derived factor-1 α , suppress amyloid β -induced neurotoxicity. **Toxicol Appl Pharmacol**. [Epub ahead of print] PMID: 21704645, 2011.
44. Rogers JT, Rusiana I, Trotter J, Zhao L, Donaldson E, Pak DT, Babus LW, Peters M, Banko JL, Chavis P, Rebeck GW, Hoe HS, Weeber EJ. Reelin supplementation enhances cognitive ability, synaptic plasticity, and dendritic spine density. **Learn Mem**. 18(9):558-64, 2011. Qin M, Xia Z, Huang T, Smith CB. Effects of chronic immobilization stress on anxiety-like behavior and basolateral amygdala morphology in Fmr1 knockout mice. **Neuroscience**. 194:282-90, 2011.
45. Rogers JT, Rusiana I, Trotter J, Zhao L, Donaldson E, Pak DT, Babus LW, Peters M, Banko JL, Chavis P, Rebeck GW, Hoe HS, Weeber EJ. Reelin supplementation enhances cognitive ability, synaptic plasticity, and dendritic spine density. **Learn Mem**. 18(9):558-64, 2011.
46. Schofield CM, Hsu R, Barker AJ, Gertz CC, Blelloch R, and Ullian EM. Monoallelic deletion of the microRNA biogenesis gene Dgcr8 produces deficits in the development of excitatory synaptic transmission in the prefrontal cortex. **Neural Dev**. 6:11, 2011.
47. Speed HE, Blaiss CA, Kim A, Haws ME, Melvin NR, Jennings M, Eisch AJ, Powell CM. Delayed reduction of hippocampal synaptic transmission and spines following exposure to repeated, subclinical doses of organophosphorus pesticide in adult mice. **Toxicol Sci**. [Epub ahead of print] PMID: 21948870, 2011.
48. Stranahan AM. Physiological variability in brain-derived neurotrophic factor expression predicts dendritic spine density in the mouse dentate gyrus. **Neurosci. Lett**. 495(1):60-2, 2011.
49. Takashima N, Odaka YS, Sakoori K, Akagi T, Hashikawa T, Morimura N, Yamada K, Aruga J. Impaired cognitive function and altered hippocampal synapse morphology in mice lacking Lrrtm1, a gene associated with schizophrenia. **PLoS One**. 6(7):e22716, 2011.

50. Tan AM, Chang YW, Zhao P, Hains BC, Waxman SG. Rac1-regulated dendritic spine remodeling contributes to neuropathic pain after peripheral nerve injury. **Exp Neurol.** 232(2):222-33, 2011.
51. Tan LA, Chawaf AA, Vaccarino FJ, Boutros PC, and Lovejoy DA. Teneurin C-terminal associated peptide (TCAP)-1 modulates dendritic morphology in hippocampal neurons and decreases anxiety-like behaviors in rats. **Physiol. & Behav.** 104(2):199-204, 2011.
52. Tao J, Wu H, Lin Q, Wei W, Lu X-H, Cantle JP, Ao Y, Olsen RW, Yang XW, Mody I, Sofroniew MV, and Sun YE. Deletion of astroglial dicer causes non-cell-autonomous neuronal dysfunction and degeneration. **J. Neurosci.** 31(22):8306-19, 2011.
53. Tereshchenko Y, Morellini F, Dityatev A, Schachner M, Irintchev A. Neural cell adhesion molecule ablation in mice causes hippocampal dysplasia and loss of septal cholinergic neurons. **J. Comp Neurol.** 519(12):2475-92, 2011.
54. Todorov B, Kros L, Shyti R, Plak P, Haasdijk ED, Raike RS, Frants RR, Hess EJ, Hoebeek FE, De Zeeuw CI, van den Maagdenberg AM. Purkinje Cell-Specific Ablation of Ca(V)2.1 Channels is Sufficient to Cause Cerebellar Ataxia in Mice. **Cerebellum.** PMID: 21870131, 2011.
55. Uchida S, Hara K, Kobayashi A, Fujimoto M, Otsuki K, Yamagata H, Hobara T, Abe N, Higuchi F, Shibata T, Hasegawa S, Kida S, Nakai A, Watanabe Y. Impaired hippocampal spinogenesis and neurogenesis and altered affective behavior in mice lacking heat shock factor 1. **Proc Natl Acad Sci U S A.** 108(4):1681-6, 2011.
56. Wang L, Budolfson K, and Wang F. Pik3c3 deletion in pyramidal neurons results in loss of synapses, extensive gliosis and progressive neurodegeneration. **Neurosci.** 172:427-42, 2011.
57. Workman JL, Manny N, Walton JC, Nelson RJ. Short day lengths alter stress and depressive-like responses, and hippocampal morphology in Siberian hamsters. **Horm Behav.** [Epub ahead of print] PMID: 21851822, 2011.
58. Xie Z, Cahill ME, Radulovic J, Wang J, Campbell SL, Miller CA, Sweatt JD, and Penzes P. Hippocampal phenotypes in kalirin-deficient mice. **Mol. Cell Neurosci.** 46(1):45-54, 2011.
59. Xu P, Das M, Reilly J, and Davis RJ. JNK regulates FoxO-dependent autophagy in neurons. **Gene Dev.** 25:310-22, 2011.
60. Yau SY, Lau BW, Tong JB, Wong R, Ching YP, Qiu G, Tang SW, Lee TM, So KF. Hippocampal neurogenesis and dendritic plasticity support running-improved spatial learning and depression-like behaviour in stressed rats. **PLoS One.** 6(9):e24263, 2011.
61. Zhang L, Yokoi F, Jin Y-H, DeAndrade MP, Hashimoto K, Standaert DG, and Li Y. Altered Dendritic Morphology of Purkinje cells in *Dyt1* ΔGAG Knock-In and Purkinje Cell-Specific *Dyt1* Conditional Knockout Mice. **PLoS One.** 6(3):e18357, 2011.
62. Zohar J, Yahalom H, Kozlovsky N, Cwikel-Hamzany S, Matar MA, Kaplan Z, Yehuda R, Cohen H. High dose hydrocortisone immediately

- after trauma may alter the trajectory of PTSD: Interplay between clinical and animal studies. **Eur Neuropsychopharmacol.** 21(11):796-809, 2011.
63. Abazyan B, Nomura J, Kannan G, Ishizuka K, Tamashiro KL, Nucifora F, Pogorelov V, Ladenheim B, Yang C, Krasnova IN, Cadet JL, Pardo C, Mori S, Kamiya A, Vogel MW, Sawa A, Ross CA, and Pletnikov MV. Prenatal interaction of mutant DISC1 and immune activation produces adult psychopathology. **Biol. Psychiat.** 68(12):1172-81, 2010.
 64. Ampuero E, Rubio FJ, Falcon R, Sandoval M, Diaz-Veliz G, Gonzalez RE, Earle N, Dagnino-Subiabre A, Aboitiz F, Orrego F, and Wyneken U. Chronic fluoxetine treatment induces structural plasticity and selective changes in glutamate receptor subunits in the rat cerebral cortex. **Neurosci.** 169(1):98-108, 2010.
 65. Ballester-Rosado CJ, Albright MJ, Wu C-S, Liao C-C, Zhu J, Xu J, Lee L-J, and Lu H-C. mGluR5 in cortical excitatory neurons exerts both cell-autonomous and -nonautonomous influences on cortical somatosensory circuit formation. **J. Neurosci.** 30(50):16896-909, 2010.
 66. Bock AS, Olavarria JF, Leigland LA, Taber EN, Jespersen SN, and Krokne CD. Diffusion tensor imaging detects early cerebral cortex abnormalities in neuronal architecture induced by bilateral neonatal enucleation: an experimental model in the ferret. **Front. Syst. Neurosci.** 4:149, 2010.
 67. Bose M, Muñoz-Illanco P, Roychowdhury S, Nichols JA, Jakkamsetti V, Porter B, Byrapureddy R, Salgado H, Kilgard MP, Aboitiz F, Dagnino-Subiabre A, and Atzori M. Effect of the environment on the dendritic morphology of the rat auditory cortex. **Synapse.** 64(2):97-110, 2010.
 68. Crupi R, Mazzon E, Marino A, La Spada G, Bramanti P, Cuzzocrea S, and Spina E. Melatonin treatment mimics the antidepressant action in chronic corticosterone-treated mice. **J. Pineal Res.** 49(2):123-9, 2010.
 69. Cubelos B, Sebastián-Serrano A, Beccari L, Calcagnotto ME, Cisneros E, Kim S, Dopazo A, Alvarez-Dolado M, Redondo JM, Bovolenta P, Walsh CA, and Nieto M. Cux1 and Cux2 regulate dendritic branching, spine morphology, and synapses of the upper layer neurons of the cortex. **Neuron.** 66(4):523-35, 2010.
 70. Deng JV, Rodriguiz RM, Hutchinson AN, Kim IH, Wetsel WC, and West AE. MeCP2 in the nucleus accumbens contributes to neural and behavioral responses to psychostimulants. **Nature Neurosci.** 13(9):1128-36, 2010.
 71. Dziras K, Coque L, Sidor MM, Kumar S, Dancy EA, Takahashi JS, McClung CA, and Nicoletis MAL. Lithium ameliorates nucleus accumbens phase-signaling dysfunction in a genetic mouse model of mania. **J. Neurosci.** 30(48):16314-23, 2010.
 72. Eiland L, and McEwen BS. Early life stress followed by subsequent adult chronic stress potentiates anxiety and blunts hippocampal structural remodeling. **Hippocampus.** doi: 10.1002/hipo.20862, 2010.
 73. Elhardt M, Martinez L, and Tejada-Simon MV. Neurochemical, behavioral and architectural changes after chronic inactivation of NMDA receptors in mice. **Neurosci. Lett.** 468(2):166-71, 2010.
 74. Fumagalli F, Pasini M, Sartorius A, Scherer R, Racagni G, Riva MA, Gass

- P. Repeated electroconvulsive shock (ECS) alters the phosphorylation of glutamate receptor subunits in the rat hippocampus. **Int J Neuropsychopharmacol.** 13(9):1255-60, 2010.
75. Hillen H, Barghorn S, Striebing A, Labkovsky B, Müller R, Nimmrich V, Nolte MW, Perez-Cruz C, van der Auwera I, van Leuven F, van Gaalen M, Beshpalov AY, Schoemaker H, Sullivan JP, and Ebert U. Generation and therapeutic efficacy of highly oligomer-specific beta-amyloid antibodies. **J. Neurosci.** 30(31):10369-79, 2010.
 76. Jentarra GM, Olfers SL, Rice SG, Srivastava N, Homanics GE, Blue M, Naidu S, and Narayanan V. Abnormalities of cell packing density and dendritic complexity in the MeCP2 A140V mouse model of Rett syndrome/X-linked mental retardation. **BMC Neurosci.** 11:19, 2010.
 77. Kelly EA, Tremblay M-E, McCasland JS, and Majewska AK. Postsynaptic deregulation in GAP-43 heterozygous mouse barrel cortex. **Cereb. Cortex.** 20(7):1696-707, 2010.
 78. Kim Y-S, Harry GJ, Kang HS, Goulding D, Wine RN, Kissling GE, Liao G, and Jetten AM. Altered cerebellar development in nuclear receptor TAK1/TR4 null mice is associated with deficits in GLAST glia, alterations in social behavior, motor learning, startle reactivity, and microglia. **Cerebellum.** 9(3):310-23, 2010.
 79. Kojima N, Hanamura K, Yamazaki H, Ikeda T, Itoharu S, and Shirao T. Genetic disruption of the alternative splicing of drebrin gene impairs context-dependent fear learning in adulthood. **Neurosci.** 165(1):138-50, 2010.
 80. Kumada T, Komuro Y, Li Y, Hu T, Wang Z, Littner Y, and Komuro H. Inhibition of cerebellar granule cell turning by alcohol. **Neurosci.** 170(4):1328-44, 2010.
 81. Langwieser N, Christel CJ, Kleppisch T, Hofmann F, Wotjak CT, and Moosmang S. Homeostatic switch in hebbian plasticity and fear learning after sustained loss of Cav1.2 calcium channels. **J. Neurosci.** 30(25):8367-75, 2010.
 82. Lee KJ, Moussa CE, Lee Y, Sung Y, Howell BW, Turner RS, Pak DT, and Hoe HS. Beta amyloid-independent role of amyloid precursor protein in generation and maintenance of dendritic spines. **Neurosci.** 169(1):344-56, 2010.
 83. Leuner B, and Gould E. Dendritic growth in medial prefrontal cortex and cognitive flexibility are enhanced during the postpartum period. **J. Neurosci.** 30(40):13499-503, 2010.
 84. Liu Q, Trotter J, Zhang J, Peters MM, Cheng H, Bao J, Han X, Weeber EJ, and Bu G. Neuronal LRP1 knockout in adult mice leads to impaired brain lipid metabolism and progressive, age-dependent synapse loss and neurodegeneration. **J. Neurosci.** 30(50):17068-78, 2010.
 85. McLaughlin KJ, Wilson JO, Harman J, Wright RL, Wiczorek L, Gomez J, Korol DL, and Conrad CD. Chronic 17 beta-estradiol or cholesterol prevents stress-induced hippocampal CA3 dendritic retraction in ovariectomized female rats: possible correspondence between CA1 spine properties and spatial acquisition. **Hippocampus.** 20(6):768-86, 2010.

86. McNair K, Spike R, Guilding C, Prendergast GC, Stone TW, Cobb SR, and Morris BJ. A role for RhoB in synaptic plasticity and the regulation of neuronal morphology. **J. Neurosci.** 30(9):3508-17, 2010.
87. Michaelson K, Zagrebelsky M, Berndt-Huch J, Polack M, Buschler A, Sendtner M, and Korte M. Neurotrophin receptors TrkB.T1 and p75NTR cooperate in modulating both functional and structural plasticity in mature hippocampal neurons. **Eur. J. Neurosci.** 32(11):1854-65, 2010.
88. Milatovic D, Montine TJ, Zaja-Milatovic S, Madison JL, Bowman AB, and Aschner M. Morphometric analysis in neurodegenerative disorders. **Curr. Protoc. Toxicol.** 46: Chapter 12, Unit 12.16.1-14, 2010.
89. Mokalled MH, Johnson A, Kim Y, Oh J, and Olson EN. Myocardin-related transcription factors regulate the Cdk5/Pctaire1 kinase cascade to control neurite outgrowth, neuronal migration and brain development. **Development.** 137(14):2365-74, 2010.
90. Nelson R. Photoperiodic regulation of affective responses and hippocampal cell morphology in Siberian hamsters. **Edward F Hayes Graduate Forum.** <http://hdl.handle.net/1811/45523>, 2010.
91. Okerlund ND, Kivimäe S, Tong CK, Peng IF, Ullian EM, and Cheyette BN. Dact1 is a postsynaptic protein required for dendrite, spine, and excitatory synapse development in the mouse forebrain. **J. Neurosci.** 30(12):4362-8, 2010.
92. Quevedo C, Sauzeau V, Menacho-Márquez M, Castro-Castro A, and Bustelo XR. Vav3-deficient mice exhibit a transient delay in cerebellar development. **Mol. Biol. Cell.** 21(6):1125-39, 2010.
93. Rebeck GW, Hoe HS, and Moussa CE. Beta-amyloid1-42 gene transfer model exhibits intraneuronal amyloid, gliosis, tau phosphorylation, and neuronal loss. **J. Biol. Chem.** 285(10):7440-6, 2010.
94. Salas-Ramirez KY, Frankfurt M, Alexander A, Luine VN, and Friedman E. Prenatal cocaine exposure increases anxiety, impairs cognitive function and increases dendritic spine density in adult rats: influence of sex. **Neurosci.** 169(3):1287-95, 2010.
95. Shmelknov SV, Hormigo A, Jing D, Proenca CC, Bath KG, Milde T, Shmelknov E, Kushner JS, Baljevic M, Dincheva I, Murphy AJ, Valenzuela DM, Gale NW, Yancopoulos GD, Ninan I, Lee FS, and Raffi S. Slitrk5 deficiency impairs corticostriatal circuitry and leads to obsessive-compulsive-like behaviors in mice. **Nature Med.** 16:598-602, 2010.
96. Sun M, Gewirtz JC, Bofenkamp L, Wickham RJ, Ge H, and O'Connor MB. Canonical TGF- β signaling is required for the balance of excitatory/inhibitory transmission within the hippocampus and prepulse inhibition of acoustic startle. **J. Neurosci.** 30(17):6025-35, 2010.
97. Wenzel HJ, Hunsaker MR, Greco CM, Willemsen R, and Berman RF. Ubiquitin-positive intranuclear inclusions in neuronal and glial cells in a mouse model of the fragile X premutation. **Brain Res.** 1318:155-66, 2010.
98. Xie Y, Hayden MR, and Xu B. BDNF overexpression in the forebrain rescues Huntington's disease phenotypes in YAC128 mice. **J. Neurosci.** 30(44):14708-18, 2010.

99. Xie Z, Cahill ME, and Penzes P. Kalirin loss results in cortical morphological alterations. **Mol. Cell Neurosci.** 43(1):81-9, 2010.
100. Akashi K, Kakizaki T, Kamiya H, Fukaya M, Yamasaki M, Abe M, Natsume R, Watanabe M, and Sakimura K. NMDA receptor GluN2B (GluR epsilon 2/NR2B) subunit is crucial for channel function, postsynaptic macromolecular organization, and actin cytoskeleton at hippocampal CA3 synapses. **J. Neurosci.** 29(35):10869-82, 2009.
101. Barros CS, Calabrese B, Chamero P, Roberts AJ, Korzus E, Lloyd K, Stowers L, Mayford M, Halpain S, and Muller U. Impaired maturation of dendritic spines without disorganization of cortical cell layers in mice lacking NRG1/ErbB signaling in the central nervous system. **Proc. Natl. Acad. Sci. USA.** 106(11):4507-12, 2009.
102. Bonett RM, Hu F, Bagamasbad P, and Denver RJ. Stressor and glucocorticoid-dependent induction of the immediate early gene *kruppel-like factor 9*: implications for neural development and plasticity. **Endocrinology.** 150(4):1757-65, 2009.
103. Cahill ME, Xie Z, Day M, Barbolina MV, Miller CA, Weiss C, Radulovic J, Sweatt JD, Disterhoft JF, Surmeier DJ, and Penzes P. Kalirin regulates cortical spine morphogenesis and disease-related behavioral phenotypes. **Proc. Natl. Acad. Sci. USA.** 106(31):13058-63, 2009.
104. Carim-Todd L, Bath KG, Fulgenzi G, Yanpallewar S, Jing D, Barrick CA, Becker J, Buckley H, Dorsey SG, Lee FS, and Tessarollo L. Endogenous truncated TrkB.T1 receptor regulates neuronal complexity and TrkB kinase receptor function in vivo. **J. Neurosci.** 29(3):678-85, 2009.
105. Chen CC, Abrams S, Pinhas A, and Brumberg JC. Morphological heterogeneity of layer VI neurons in mouse barrel cortex. **J. Comp. Neurol.** 512(6):726-46, 2009.
106. Chen ZL, Haegeli V, Yu H, and Strickland S. Cortical deficiency of laminin gamma1 impairs the AKT/GSK-3beta signaling pathway and leads to defects in neurite outgrowth and neuronal migration. **Dev. Biol.** 327(1):158-68, 2009.
107. Dalla C, Whetstone AS, Hodes GE, and Shors TJ. Stressful experience has opposite effects on dendritic spines in the hippocampus of cycling versus masculinized females. **Neurosci. Lett.** 449(1):52-6, 2009.
108. Dumanis SB, Tesoriero JA, Babus LW, Nguyen MT, Trotter JH, Ladu MJ, Weeber EJ, Turner RS, Xu B, Rebeck GW, and Hoe HS. ApoE4 decreases spine density and dendritic complexity in cortical neurons in vivo. **J. Neurosci.** 29(48):15317-22, 2009.
109. Enard W, Gehre S, Hammerschmidt K, Hölter SM, Blass T, Somel M, Brückner MK, Schreiweis C, Winter C, Sohr R, Becker L, Wiebe V, Nickel B, Giger T, Müller U, Groszer M, Adler T, Aguilar A, Bolle I, Calzada-Wack J, Dalke C, Ehrhardt N, Favor J, Fuchs H, Gailus-Durner V, Hans W, Hölzlwimmer G, Javaheri A, Kalaydjiev S, Kallnik M, Kling E, Kunder S, Mossbrugger I, Naton B, Racz I, Rathkolb B, Rozman J, Schrewe A, Busch DH, Graw J, Ivandic B, Klingenspor M, Klopstock T, Ollert M, Quintanilla-Martinez L, Schulz H, Wolf E, Wurst W, Zimmer A,

- Fisher SE, Morgenstern R, Arendt T, de Angelis MH, Fischer J, Schwarz J, and Pääbo S. A humanized version of Foxp2 affects cortico-basal ganglia circuits in mice. **Cell**. 137(5):961-71, 2009.
110. Frankfurt M, Wang HY, Marmolejo N, Bakshi K, and Friedman E. Prenatal cocaine increases dendritic spine density in cortical and subcortical brain regions of the rat. **Dev. Neurosci**. 31(1-2):71-5, 2009.
 111. Goshen I, Avital A, Kreisel T, Licht T, Segal M, and Yirmiya R. Environmental enrichment restores memory functioning in mice with impaired IL-1 signaling via reinstatement of long-term potentiation and spine size enlargement. **J. Neurosci**. 29(11):3395-403, 2009.
 112. Griffin GD, and Flanagan-Cato LM. Sex differences in the dendritic arbor of hypothalamic ventromedial nucleus neurons. **Physiol. Behav.** 97(2):151-6, 2009.
 113. Hains AB, Vu MA, Maciejewski PK, van Dyck CH, Gotttron M, and Arnsten AF. Inhibition of protein kinase C signaling protects prefrontal cortex dendritic spines and cognition from the effects of chronic stress. **Proc. Natl. Acad. Sci. USA**. 106(42):17957-62, 2009.
 114. Hoe HS, Lee KJ, Carney RS, Lee J, Markova A, Lee JY, Howell BW, Hyman BT, Pak DT, Bu G, and Rebeck GW. Interaction of reelin with amyloid precursor protein promotes neurite outgrowth. **J. Neurosci**. 29(23):7459-73, 2009.
 115. Hoskison MM, Moore AN, Hu B, Orsi S, Kobori N, and Dash PK. Persistent working memory dysfunction following traumatic brain injury: evidence for a time-dependent mechanism. **Neurosci**. 159(2):483-91, 2009.
 116. Jakovcevski I, Siering J, Hargus G, Karl N, Hoelters L, Djogo N, Yin S, Zecevic N, Schachner M, and Irintchev A. Close homologue of adhesion molecule L1 promotes survival of Purkinje and granule cells and granule cell migration during murine cerebellar development. **J. Comp. Neurol**. 513(5):496-510, 2009.
 117. Johnson SA, Wang JF, Sun X, McEwen BS, Chattarji S, and Young LT. Lithium treatment prevents stress-induced dendritic remodeling in the rodent amygdala. **Neurosci**. 163(1):34-9, 2009.
 118. Karlén A, Karlsson TE, Mattsson A, Lundströmer K, Codeluppi S, Pham TM, Bäckman CM, Ogren SO, Aberg E, Hoffman AF, Sherling MA, Lupica CR, Hoffer BJ, Spenger C, Josephson A, Brené S, Olson L. Nogo receptor 1 regulates formation of lasting memories. **Proc Natl Acad Sci U S A**. 106(48):20476-81, 2009.
 119. Kerjan G, Koizumi H, Han EB, Dubé CM, Djakovic SN, Patrick GN, Baram TZ, Heinemann SF, and Gleeson JG. Mice lacking doublecortin and doublecortin-like kinase 2 display altered hippocampal neuronal maturation and spontaneous seizures. **Proc. Natl. Acad. Sci. USA**. 106(16):6766-71, 2009.
 120. Kumada T, Jiang Y, Kawanami A, Cameron DB, and Komuro H. Autonomous turning of cerebellar granule cells in vitro by intrinsic programs. **Dev. Biol**. 326(1):237-49, 2009.
 121. Labelle DR, Cox JM, Dunn-Meynell AA, Levin BE, and Flanagan-Cato

- LM. Genetic and dietary effects on dendrites in the rat hypothalamic ventromedial nucleus. **Physiol. Behav.** 98(4):511-6, 2009.
122. Melendez-Ferro M, Perez-Costas E, and Roberts RC. A new use for long-term frozen brain tissue: golgi impregnation. **J. Neurosci. Methods.** 176(2):72-7, 2009.
 123. Ng D, Pitcher GM, Szilard RK, Sertié A, Kanisek M, Clapcote SJ, Lipina T, Kalia LV, Joo D, McKerlie C, Cortez M, Roder JC, Salter MW, and McInnes RR. Neto1 is a novel CUB-domain NMDA receptor-interacting protein required for synaptic plasticity and learning. **PLoS Biol.** 7(2):e41, 2009.
 124. Nithianantharajah J, Barkus C, Vijiaratnam N, Clement O, and Hannan AJ. Modeling brain reserve: experience-dependent neuronal plasticity in healthy and Huntington's disease transgenic mice. **Am. J. Geriatr. Psychiat.** 17(3):196-209, 2009.
 125. Rajasekaran K, Sun C, and Bertram EH. Altered pharmacology and GABA-A receptor subunit expression in dorsal midline thalamic neurons in limbic epilepsy. **Neurobiol. Dis.** 33(1):119-32, 2009.
 126. Schaefer A, Sampath SC, Intrator A, Min A, Gertler TS, Surmeier DJ, Tarakhovsky A, and Greengard P. Control of cognition and adaptive behavior by the GLP/G9a epigenetic suppressor complex. **Neuron.** 64(5):678-91, 2009.
 127. Scobie KN, Hall BJ, Wilke SA, Klemenhausen KC, Fujii-Kuriyama Y, Ghosh A, Hen R, and Sahay A. Krüppel-like factor 9 is necessary for late-phase neuronal maturation in the developing dentate gyrus and during adult hippocampal neurogenesis. **J. Neurosci.** 29(31):9875-87, 2009.
 128. Shi Q, Hu X, Prior M, and Yan R. The occurrence of aging-dependent reticulum 3 immunoreactive dystrophic neurites decreases cognitive function. **J. Neurosci.** 29(16):5108-15, 2009.
 129. Stranahan AM, Lee K, Martin B, Maudsley S, Golden E, Cutler RG, and Mattson MP. Voluntary exercise and caloric restriction enhance hippocampal dendritic spine density and BDNF levels in diabetic mice. **Hippocampus.** 19(10):951-61, 2009.
 130. Tan AM, Choi JS, Waxman SG, and Hains BC. Dendritic spine remodeling after spinal cord injury alters neuronal signal processing. **J. Neurophysiol.** 102(4):2396-409, 2009.
 131. Tran TS, Rubio ME, Clem RL, Johnson D, Case L, Tessier-Lavigne M, Huganir RL, Ginty DD, and Kolodkin AL. Secreted semaphorins control spine distribution and morphogenesis in the postnatal CNS. **Nature.** 462(7276):1065-9, 2009.
 132. Tropea D, Giacometti E, Wilson NR, Beard C, McCurry C, Fu DD, Flannery R, Jaenisch R, and Sur M. Partial reversal of Rett Syndrome-like symptoms in MeCP2 mutant mice. **Proc. Natl. Acad. Sci. USA.** 106(6):2029-34, 2009.
 133. Ullner PM, Di Nardo A, Goldman JE, Schobel S, Yang H, Engelstad K, Wang D, Sahin M, and De Vivo DC. Murine Glut-1 transporter haploinsufficiency: postnatal deceleration of brain weight and reactive astrocyto-

- sis. **Neurobiol. Dis.** 36(1):60-9, 2009.
134. Wang HD, Stanwood GD, Grandy DK, and Deutch AY. Dystrophic dendrites in prefrontal cortical pyramidal cells of dopamine D1 and D2 but not D4 receptor knockout mice. **Brain Res.** 1300:58-64, 2009.
 135. Workman JL, Bowers SL, and Nelson RJ. Enrichment and photoperiod interact to affect spatial learning and hippocampal dendritic morphology in white-footed mice (*Peromyscus leucopus*). **Eur. J. Neurosci.** 29(1):161-70, 2009.
 136. Xu Y, Lin D, Li S, Li G, Shyamala SG, Barish PA, Vernon MM, Pan J, and Ogle WO. Curcumin reverses impaired cognition and neuronal plasticity induced by chronic stress. **Neuropharmacology.** 57(4):463-71, 2009.
 137. Yu H, Wang Y, Pattwell S, Jing D, Liu T, Zhang Y, Bath KG, Lee FS, and Chen ZY. Variant BDNF Val66Met polymorphism affects extinction of conditioned aversive memory. **J. Neurosci.** 29(13):4056-64, 2009.
 138. An JJ, Gharami K, Liao GY, Woo NH, Lau AG, Vanevski F, Torre ER, Jones KR, Feng Y, Lu B, and Xu B. Distinct role of long 3' UTR BDNF mRNA in spine morphology and synaptic plasticity in hippocampal neurons. **Cell.** 134(1):175-87, 2008.
 139. Brun VH, Leutgeb S, Wu HQ, Schwarcz R, Witter MP, Moser EI, and Moser MB. Impaired spatial representation in CA1 after lesion of direct input from entorhinal cortex. **Neuron.** 57(2):290-302, 2008.
 140. Chen Q, Xiong X, Lee TH, Liu Y, Sun QA, Wetsel W, and Zhang X. Adeno-associated virus-mediated ILK gene silencing in the rat NAc core. **J. Neurosci. Methods.** 173(2):208-14, 2008.
 141. Chen Q, Xiong X, Lee TH, Liu Y, Wetsel WC, and Zhang X. Neural plasticity and addiction: integrin-linked kinase and cocaine behavioral sensitization. **J. Neurochem.** 107(3):679-89, 2008.
 142. Davis TH, Cuellar TL, Koch SM, Barker AJ, Harfe BD, McManus MT, and Ullian EM. Conditional loss of Dicer disrupts cellular and tissue morphogenesis in the cortex and hippocampus. **J. Neurosci.** 28(17):4322-30, 2008.
 143. Dindot SV, Antalffy BA, Bhattacharjee MB, and Beaudet AL. The Angelman syndrome ubiquitin ligase localizes to the synapse and nucleus, and maternal deficiency results in abnormal dendritic spine morphology. **Hum. Mol. Genet.** 17(1):111-8, 2008.
 144. Fraser MM, Bayazitov IT, Zakharenko SS, and Baker SJ. Phosphatase and tensin homolog, deleted on chromosome 10 deficiency in brain causes defects in synaptic structure, transmission and plasticity, and myelination abnormalities. **Neurosci.** 151(2):476-88, 2008.
 145. Griffin GD, and Flanagan-Cato LM. Estradiol and progesterone differentially regulate the dendritic arbor of neurons in the hypothalamic ventromedial nucleus of the female rat (*Rattus norvegicus*). **J. Comp. Neurol.** 510(6):631-40, 2008.
 146. Kosodo Y, Toida K, Dubreuil V, Alexandre P, Schenk J, Kiyokage E, Attardo A, Mora-Bermúdez F, Arii T, Clarke JD, and Huttner WB. Cytokinesis of neuroepithelial cells can divide their basal process before ana-

- phase. **EMBO J.** 27(23):3151-63, 2008.
147. Kurz JE, Moore BJ, Henderson SC, Campbell JN, and Churn SB. A cellular mechanism for dendritic spine loss in the pilocarpine model of status epilepticus. **Epilepsia.** 49(10):1696-710, 2008.
 148. Leão RN, Leão RM, da Costa LF, Rock Levinson S, and Walmsley B. A novel role for MNTB neuron dendrites in regulating action potential amplitude and cell excitability during repetitive firing. **Eur. J. Neurosci.** 27(12):3095-108, 2008.
 149. Li CQ, Liu D, Huang L, Wang H, Zhang JY, and Luo XG. Cytosine arabinoside treatment impairs the remote spatial memory function and induces dendritic retraction in the anterior cingulate cortex of rats. **Brain Res. Bull.** 77(5):237-40, 2008.
 150. Li X, Zhang P, Yang Y, Xiong Y, Qi Y, and Hu H. Differentiation and developmental origin of cerebellar granule neuron ectopia in protein O-mannose UDP-N-acetylglucosaminyl transferase 1 knockout mice. **Neurosci.** 152(2):391-406, 2008.
 151. Liu QA, and Shio H. Mitochondrial morphogenesis, dendrite development, and synapse formation in cerebellum require both Bcl-w and the glutamate receptor delta2. **PLoS Genet.** 4(6):e1000097, 2008.
 152. Mishra A, Knerr B, Paixão S, Kramer ER, and Klein R. The protein dendrite arborization and synapse maturation 1 (Dasm-1) is dispensable for dendrite arborization. **Mol. Cell Biol.** 28(8):2782-91, 2008.
 153. Pandey SC, Zhang H, Ugale R, Prakash A, Xu T, and Misra K. Effector immediate-early gene arc in the amygdala plays a critical role in alcoholism. **J. Neurosci.** 28(10):2589-600, 2008.
 154. Papadeas ST, Halloran C, McCown TJ, Breese GR, and Blake BL. Changes in apical dendritic structure correlate with sustained ERK1/2 phosphorylation in medial prefrontal cortex of a rat model of dopamine D1 receptor agonist sensitization. **J. Comp. Neurol.** 511(2):271-85, 2008.
 155. Quach TT, Massicotte G, Belin MF, Honnorat J, Glasper ER, Devries AC, Jakeman LB, Baudry M, Duchemin AM, and Kolattukudy PE. CRMP3 is required for hippocampal CA1 dendritic organization and plasticity. **FASEB J.** 22(2):401-9, 2008.
 156. Samuels IS, Karlo JC, Faruzzi AN, Pickering K, Herrup K, Sweatt JD, Saitta SC, and Landreth GE. Deletion of ERK2 mitogen-activated protein kinase identifies its key roles in cortical neurogenesis and cognitive function. **J. Neurosci.** 28(27):6983-95, 2008.
 157. Savonenko AV, Melnikova T, Laird FM, Stewart KA, Price DL, and Wong PC. Alteration of BACE1-dependent NRG1/ErbB4 signaling and schizophrenia-like phenotypes in BACE1-null mice. **Proc. Natl. Acad. Sci. USA.** 105(14):5585-90, 2008.
 158. Tan AM, Stambouliau S, Chang YW, Zhao P, Hains AB, Waxman SG, and Hains BC. Neuropathic pain memory is maintained by Rac1-regulated dendritic spine remodeling after spinal cord injury. **J. Neurosci.** 28(49):13173-83, 2008.
 159. Yamasaki N, Maekawa M, Kobayashi K, Kajii Y, Maeda J, Soma M,

- Takao K, Tanda K, Ohira K, Toyama K, Kanzaki K, Fukunaga K, Sudo Y, Ichinose H, Ikeda M, Iwata N, Ozaki N, Suzuki H, Higuchi M, Suhara T, Yuasa S, and Miyakawa T. Alpha-CaMKII deficiency causes immature dentate gyrus, a novel candidate endophenotype of psychiatric disorders. **Mol. Brain.** 1(1):6, 2008.
160. Ye H, Tan YL, Ponniah S, Takeda Y, Wang SQ, Schachner M, Watanabe K, Pallen CJ, and Xiao ZC. Neural recognition molecules CHL1 and NB-3 regulate apical dendrite orientation in the neocortex via PTP alpha. **EMBO J.** 27(1):188-200, 2008.
 161. Zaja-Milatovic S, Gupta RC, Aschner M, Montine TJ, and Milatovic D. Pharmacologic suppression of oxidative damage and dendritic degeneration following kainic acid-induced excitotoxicity in mouse cerebrum. **Neurotoxicol.** 29(4):621-7, 2008.
 162. Zhang L, Hernández VS, Medina-Pizarro M, Valle-Leija P, Vega-González A, and Morales T. Maternal hyperthyroidism in rats impairs stress coping of adult offspring. **J. Neurosci. Res.** 86(6):1306-15, 2008.
 163. Ampuero E, Dagnino-Subiabre A, Sandoval R, Zepeda-carreno R, Sandoval S, Viedma A, Aboitiz F, Orrego F, and Wyneken U. Status epilepticus induces region-specific changes in dendritic spines, dendritic length and TrkB protein content of rat brain cortex. **Brain Res.** 1150:225-38, 2007.
 164. Conrad CD, McLaughlin KJ, Harman JS, Foltz C, Wiczorek L, Lightner E, and Wright RL. Chronic glucocorticoids increase hippocampal vulnerability to neurotoxicity under conditions that produce CA3 dendritic retraction but fail to impair spatial recognition memory. **J. Neurosci.** 27:8278-85, 2007.
 165. Cornejo BJ, Mesches MH, Coultrap S, Browning MD, and Benke TA. A single episode of neonatal seizures permanently alters glutamatergic synapses. **Ann. Neurol.** 61:411-26, 2007.
 166. Fünfschilling U, Saher G, Xiao L, Möbius W, and Nave A. Survival of adult neurons lacking cholesterol synthesis in vivo. **BMC Neurosci.** 8:1, 2007.
 167. Gunnarsen JM, Kim MH, Fuller SJ, De Silva M, Britto JM, Hammond VE, Davies PJ, Petrou S, Faber ES, Sah P, and Tan SS. Sez-6 proteins affect dendritic arborization patterns and excitability of cortical pyramidal neurons. **Neuron.** 56(4):621-39, 2007.
 168. Johnson DA, Zhang J, Frase S, Wilson M, Rodriguez-Galindo C, and Dyer MA. Neuronal differentiation and synaptogenesis in retinoblastoma. **Cancer Res.** 67:2701-11, 2007.
 169. Keays DA, Tian G, Poirier K, Huang GJ, Siebold C, Cleak J, Oliver PL, Fray M, Harvey RJ, Molnar Z, Pinon MC, Dear N, Valdar W, Brown SDM, Davies KE, Rawlins JNP, Cowan NJ, Nolan P, Chelly J, and Flint J. Mutations in α -tubulin cause abnormal neuronal migration and lissencephaly in humans. **Cell.** 128:45-57, 2007.
 170. Kennedy NJ, Martin G, Ehrhardt AG, Cavanagh-Kyros J, Kuan CY, Rakic P, Flavell RA, Treisman SN, and Davis RJ. Requirement of JIP scaffold

- proteins for NMDA-mediated signal transduction. **Gene Dev.** 21:2336-46, 2007.
171. Kikkawa YS, and Pawlowski KS. Cochlear neuronal tracing for frequency mapping with DiI, NeuroVue, and Golgi methods. **Acta Oto-laryngol Suppl.** 559:19-23, 2007.
 172. Kimura T, Yamashita S, Fukuda T, Park JM, Murayama M, Mizoroki T, Yoshiike Y, Sahara N, and Takashima A. Hyperphosphorylated tau in parahippocampal cortex impairs place learning in aged mice expressing wild-type human tau. **EMBO J.** 26(24):5143-52, 2007.
 173. Li L, Yun SH, Keblesh J, Trommer BL, Xiong H, Radulovic J, and Tourtellotte WG. Egr3, a synaptic activity regulated transcription factor that is essential for learning and memory. **Mol. and Cell. Neurosci.** 35:76-88, 2007.
 174. McLaughlin KJ, Gomez JL, Baran SE, and Conrad CD. The effects of chronic stress on hippocampal morphology and function: an evaluation of chronic restraint paradigms. **Brain Res.** 1161:56-64, 2007.
 175. Nielsen JV, Nielsen FH, Ismail R, Noraberg J, and Jensen NA . Hippocampus-like corticoneurogenesis induced by two isoforms of the BTB-zinc finger gene Zbtb20 in mice. **Development.** 134(6):1133-40, 2007.
 176. Ostwald D, Karpac J, and Hochgeschwender U. Effects on hippocampus of lifelong absence of glucocorticoids in the pro-opiomelanocortin null mutant mouse reveal complex relationship between glucocorticoids and hippocampal structure and function. **J. Mol. Neurosci.** 28:291-302, 2007.
 177. Sadakata T, Kakegawa W, Mizoguchi A, Washida M, Katoh-Semba R, Shutoh F, Okamoto T, Nakashima H, Kimura K, Tanaka M, Sekine Y, Itohara S, Yuzaki M, Nagao S, and Furuichi T. Impaired cerebellar development and function in mice lacking CAPS2, a protein involved in neurotrophin release. **J. Neurosci.** 27:2472-82, 2007.
 178. Sarti F, Borgland SL, Kharazia VN, and Bonci A. Acute cocaine exposure alters spine density and long-term potentiation in the ventral tegmental area. **Eur. J. Neurosci.** 26:749-56, 2007.
 179. Shu Y, Duque A, Yu Y, Haider B, and McCormick DA. Properties of action potential initiation in neocortical pyramidal cells: evidence from whole cell axon recordings. **J. Neurophysiol.** 97:746-60, 2007.
 180. Soderling SH, Guire ES, Kaech S, White J, Zhang F, Schutz K, Langeberg LK, Banker G, Raber J, and Scott JD. A WAVE-1 and WRP signaling complex regulates spine density, synaptic plasticity, and memory. **J. Neurosci.** 27:355-65, 2007.
 181. Sun M, Thomas MJ, Herder R, Bofenkamp ML, Selleck SB, and O'Connor MB. Presynaptic contributions of chordin to hippocampal plasticity and spatial learning. **J. Neurosci.** 27:7740-50, 2007.
 182. Wallace M, Frankfurt M, Arellanos A, Inagaki T, and Luine V. Impaired recognition memory and decreased prefrontal cortex spine density in aged female rats. **Ann. N. Y. Acad. Sci.** 1097: 54-7, 2007.
 183. Watase K, Gatchel JR, Sun Y, Emamian E, Atkinson R, Richman R, Mizusawa H, Orr HT, Shaw C, and Zoghbi HY. Lithium therapy improves

- neurological function and hippocampal dendritic arborization in a spinocerebellar ataxia type 1 mouse model. **PLoS Med.** 4:836-47, 2007.
184. Adhami F, Liao GH, Morozov YM, Schloemer A, Schmithorst VJ, Lorenz JN, Dunn RS, Vorhees CV, Wills-Karp M, Degen JL, Davis RJ, Mizushima N, Rakic P, Dardzinski BJ, Holland SK, Sharp FR, and Kuan CY. Cerebral ischemia-hypoxia induces intravascular coagulation and autophagy. **Am. J. Pathol.** 169:566-83, 2006.
 185. Buss RR, Gould TW, Ma J, Vinsant S, Prevette D, Winseck A, Toops KA, Hammarback JA, Smith TL, and Oppenheim RW. Neuromuscular development in the absence of programmed cell death: phenotypic alteration of motoneurons and muscle. **J. Neurosci.** 26:13413-27, 2006.
 186. Dang MT, Yokoi FY, Yin HH, Lovinger DM, Wang Y, and Li Y. Disrupted motor learning and long-term synaptic plasticity in mice lacking NMDAR1 in the striatum. **Proc. Natl. Acad. Sci. USA.** 103:15254-9, 2006.
 187. Elia LP, Yamamoto M, Zang K, and Reichardt LF. p120 catenin regulates dendritic spine and synapse development through Rho-family GTPases and cadherins. **Neuron.** 51:43-56, 2006.
 188. Huang Z, Shimazu K, Woo NH, Zang K, Müller U, Lu B, and Reichardt LF. Distinct roles of the β 1-class integrins at the developing and the mature hippocampal excitatory synapse. **J. Neurosci.** 26:11208-19, 2006.
 189. Inan M, Lu HC, Albright MJ, She WC, and Crair MC. Barrel map development relies on protein kinase A regulatory subunit II β -mediated cAMP signaling. **J. Neurosci.** 26:4338-49, 2006.
 190. Jacobsen JS, Wu CC, Redwine JM, Comery TA, Arias R, Bowlby M, Martone R, Morrison JH, Pangalos MN, Reinhart PH, and Bloom FE. Early-onset behavioral and synaptic deficits in a mouse model of Alzheimer's disease. **Proc. Natl. Acad. Sci. USA.** 103:5161-6, 2006.
 191. Kleen JK, Sitomer MT, Killeen PR, and Conrad CD. Chronic stress impairs spatial memory and motivation for reward without disrupting motor ability and motivation to explore. **Behav. Neurosci.** 120:842-51, 2006.
 192. Luine V, Attalla S, Mohan G, Costa A, and Frankfurt M. Dietary phytoestrogens enhance spatial memory and spine density in the hippocampus and prefrontal cortex of ovariectomized rats. **Brain Res.** 1126:183-7, 2006.
 193. Maezawa I, Zaja-Milatovic S, Milatovic D, Stephen C, Sokal I, Maeda N, Montine TJ, and Montine KS. Apolipoprotein E isoform-dependent dendritic recovery of hippocampal neurons following activation of innate immunity. **J. Neuroinflamm.** 3:21, 2006.
 194. Melendez-Ferro M, Perez-Costas E, and Roberts RC. Golgi staining of long-term stored human brain tissue. **FENS Abstr.** 3:A059.20, 2006.
 195. Moretti P, Levenson JM, Battaglia F, Atkinson R, Teague R, Antalffy B, Armstrong D, Arancia O, Sweatt JD, and Zoghbi HY. Learning and memory and synaptic plasticity are impaired in a mouse model of Rett syndrome. **J. Neurosci.** 26:319-27, 2006.
 196. Poet M, Kornak U, Schweizer M, Zdebik AA, Scheel O, Hoelter S, Wurst W, Schmitt A, Fuhrmann JC, Planells-Cases R, Mole SE, Hubner CA, and

- Jentsch TJ. Lysosomal storage disease upon disruption of the neuronal chloride transport protein CIC-6. **Proc. Natl. Acad. Sci. USA.** 103:13854-9, 2006.
197. Wallace M, Luine V, Arellanos A, and Frankfurt M. Ovariectomized rats show decreased recognition memory and spine density in the hippocampus and prefrontal cortex. **Brain Res.** 1126:176-82, 2006.
 198. Watson KK, Jones TK, and Allman JM. Dendritic architecture of the von Economo neurons. **Neurosci.** 141:1107-12, 2006.
 199. Björklom B, Östman N, Hongisto V, Komarovski V, Filén JJ, Nyman TA, Kallunki T, Courtney MJ, and Coffey ET. Constitutively active cytoplasmic c-jun N-terminal kinase 1 is a dominant regulator of dendritic architecture: role of microtubule-associated protein 2 as an effector. **J. Neurosci.** 25:6350-61, 2005.
 200. Dagnino-Subiabre A, Terreros G, Carmona-Fontaine C, Zepeda R, Orellana JA, Diaz-Veliz G, Mora S, and Aboitiz F. Chronic stress impairs acoustic conditioning more than visual conditioning in rats: morphological and behavioural evidence. **Neurosci.** 135:1067-74, 2005.
 201. Grillet N, Pattyn A, Contet C, Kieffer BL, Goridis C, and JF Brunet. Generation and characterization of Rgs4 mutant mice. **Mol. and Cell. Biol.** 25:4221-8, 2005.
 202. Gu X, Li C, Wei W, Lo V, Gong S, Li SH, Iwasato T, Itoharu S, Li XJ, Mody I, Heintz N, and Yang XW. Pathological cell-cell interactions elicited by a neuropathogenic form of mutant huntingtin contribute to cortical pathogenesis in HD mice. **Neuron.** 46:433-44, 2005.
 203. Ishikura N, Clever JL, Bouzamondo-Bernstein E, Samayoa E, Prusiner SB, Huang EJ, and DeArmond SJ. Notch-1 activation and dendritic atrophy in prion disease. **Proc. Natl. Acad. Sci. USA.** 102:886-91, 2005.
 204. Laub F, Lei L, Sumiyoshi H, Kajimura D, Dragomir C, Smaldone S, Puche AC, Petros TJ, Mason C, Parada LF, and Ramirez F. Transcription factor KLF7 is important for neuronal morphogenesis in selected regions of the nervous system. **Mol. and Cell. Biol.** 25:5699-711, 2005.
 205. McLaughlin KJ, Baran SE, Wright RL, and Conrad CD. Chronic stress enhances spatial memory in ovariectomized female rats despite CA3 dendritic retraction: possible involvement of CA1 neurons. **Neurosci.** 135:1045-54, 2005.
 206. Niewmierzycka A, Mills J, St-Arnaud R, Dedhar S, and Reichardt LF. Integrin-linked kinase deletion from mouse cortex results in cortical lamination defects resembling cobblestone lissencephaly. **J. Neurosci.** 25:7022-31, 2005.
 207. Pyter LM, Reader BF, and Nelson RJ. Short photoperiods impair spatial learning and alter hippocampal dendritic morphology in adult male white-footed mice (*Peromyscus leucopus*). **J. Neurosci.** 25:4521-6, 2005.
 208. Ramanan N, Shen Y, Sarsfield S, Lemberger T, Schütz G, Linden DJ, and Ginty DD. SRF mediates activity-induced gene expression and synaptic plasticity but not neuronal viability. **Nature Neurosci.** 8:759-67, 2005.
 209. Ren-Patterson RF, Cochran LW, Holmes A, Sherrill S, Huang SJ, Tolliver

- T, Lesch KP, Lu B, and Murphy DL. Loss of brain-derived neurotrophic factor gene allele exacerbates brain monoamine deficiencies and increases stress abnormalities of serotonin transporter knockout mice. **J. Neurosci. Res.** 79:756-71, 2005.
210. Spiga S, Puddu MC, Pisano M, Diana M. Morphine withdrawal-induced morphological changes in the nucleus accumbens. **Eur J Neurosci.** 22 (9):2332-40, 2005.
 211. Amateau SK, and McCarthy MM. Induction of PGE₂ by estradiol mediates developmental masculinization of sex behavior. **Nature Neurosci.** 7:643-50, 2004.
 212. Mlatovic D, Zaja-Milatovic S, Montine KS, Horner PJ, and Montine TJ. Pharmacologic suppression of neuronal oxidative damage and dendritic degeneration following direct activation of glial innate immunity in mouse cerebrum. **J. Neurochem.** 87:1518-26, 2003.
 213. Dahl JP, Wang-Dunlop J, Gonzales C, Goad MEP, Mark RJ, and Kwak SP. Characterization of the WAVE1 knock-out mouse: implications for CNS development. **J. Neurosci.** 23:3343-52, 2003.
 214. Beggs HE, Schahin-Reed D, Zang K, Goebbels S, Nave KA, Gorski J, Jones KR, Sretavan D, and Reichardt LF. FAK deficiency in cells contributing to the basal lamina results in cortical abnormalities resembling congenital muscular dystrophies. **Neuron.** 40:501-14, 2003.

**For more references, please visit our website at
www.fdnurotech.com.**

Notes
