

## Rhod-4

*Rhod-4 is the new generation of calcium assay red dye*

### Product Information

Product name	cat.number	Packaging interest
<b>Rhod-4 AM</b>		
CQ6061	5 x 50 µg	Cell-permeable
CQ6062	10 x 50 µg	
CQ6063	20 x 50 µg	
CQ6064	1 mg	
<b>Rhod-4, sodium salt</b>		
CQ6070	10 x 50 µg	
<b>Rhod-4 NW Calcium Assay Kit, Medium Removal</b>		
CQ6080	1 plate	Performed the assay in a convenient 96-well or 384-well microtiter-plate format and easily adapted to automation.
CQ6081	10 plates	
CQ6082	100 plates	
<b>Rhod-4 NW Calcium Assay Kit, 1% FBS Growth Medium</b>		
CQ6090	1 plate	Performed the assay in a convenient 96-well or 384-well microtiter-plate format and easily adapted to automation. Higher dye concentration to avoid medium removal.
CQ6091	10 plates	
CQ6092	100 plates	
<b>Fluorescence:</b>	$\lambda_{exc./em.}$ : 530 / 555 nm	
<b>Kd (calcium)</b>	525 nM	

**Storage:** -20°C. Protect from light.

### Introduction

Calcium acts as a universal second messenger in a variety of cells. The beginning of life, the act of fertilization, is regulated by  $Ca^{2+}$ . Numerous functions of all types of cells are regulated by  $Ca^{2+}$  to a greater or lesser degree. Since the 1920s, scientists have attempted to measure  $Ca^{2+}$ , but few were successful due to limited availability of  $Ca^{2+}$  probes. The first reliable measurements of  $Ca^{2+}$  were performed by Ridgway and Ashley by injecting the photoprotein aequorin into the giant muscle fiber of the barnacle. Subsequently, in the 1980s, Tsien and colleagues produced a variety of fluorescent indicators. Among them the rhodamine-based  $Ca^{2+}$  reagents (such as Rhod-2) have been the most valuable longer wavelength dye for measuring  $Ca^{2+}$  with red fluorescence.

### Key Features

Although Rhod-2 has been the most popular red fluorescent  $Ca^{2+}$  indicator, its mitochondrial localization and high basal  $Ca^{2+}$  signal in cells have severely limited its cellular applications. In addition, the less optimal excitation of Rhod-2 at 488 nm makes it less robust to use with some instruments (such as FLIPR™) that have only 488 nm excitation light source. Our Rhod-4™ serial calcium detection reagents have been developed to address these limitations of Rhod-2.

The absorption and emission peaks of Rhod-4™ reagents are 530 nm and 555 nm, respectively. Although Rhod-4 has maximum absorption at 530 nm, its absorption at 488 nm is quite strong (see Figure 1). It is quite unique that Rhod-4™ can be well excited with an argon ion laser at 488 nm besides the longer wavelength 514 nm, 532 nm and 546 nm excitations. Rhod-4™ emits fluorescence (at wavelengths 555 nm) that increases with increasing  $Ca^{2+}$ . Rhod-4™ is determined to undergo a >200-fold increase in fluorescence upon binding  $Ca^{2+}$ . Because the range of increase in  $Ca^{2+}$  in many cells after stimulation is generally 5- to 10-fold, Rhod-4™ is an excellent probe to use with high sensitivity in this re-

[Info@fluoprobes.com](mailto:Info@fluoprobes.com)

[Technical-](#)

[support@fluoprobes.com](mailto:support@fluoprobes.com)

[Order-online@fluoprobes.com](mailto:Order-online@fluoprobes.com)

Contact your local distributor

FluoProbes®, powered by



213 Avenue J.F. Kennedy - BP 1140  
93103 Montluçon Cedex - France  
Tél. 04 70 03 88 55 - Fax 04 70 03 82 60

P.1

FT-CQ6062

gion. The  $K_d$  of Rhod-4™ is estimated to be 525 nM (22°C, pH 7.0–7.5), but this value may be significantly influenced by pH, viscosity, and binding proteins in vivo conditions


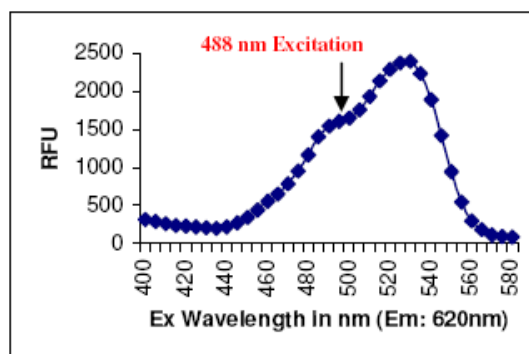


Figure 1. Excitation Spectrum of Rhod-4™, sodium salt in the presence of calcium chloride.


Besides their convenient excitation wavelengths and large fluorescence enhancement by calcium, Rhod-4™ is much brighter in cells than Rhod-2 as shown in Figure 2. More importantly Rhod-4™ is predominantly localized in cytosols unlike Rhod-2 that is mainly localized in mitochondria. In addition, Rhod-4 AM is much more readily loaded into live cells than Rhod-2 AM. Rhod-4™ reagents have a less temperature-dependent cell loading property, giving similar results either at room temperature or 37°C. This characteristic makes Rhod-4™ more robust for HTS applications than Rhod-2 AM.

Compared to Rhod-2, our Rhod-4™ calcium detection reagents have the following advantages:

- *Convenient excitation Wavelengths:* multiple excitation options @ ~490 nm, 514 nm, 532 nm and 546 nm.
- *Much Larger Assay Window:* 10 times better than Rhod-2 AM
- *Enhanced Intensity:* 4 times brighter than Rhod-2 AM
- *Faster Loading:* dye loading at room temperature
- *Versatile Packing Sizes to Meet Your Special Needs:* 1 mg; 10x50 µg; 20x50 µg; HTS packages.

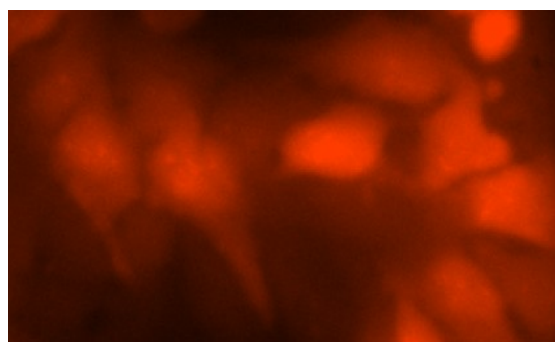

Rhod-4™ AM

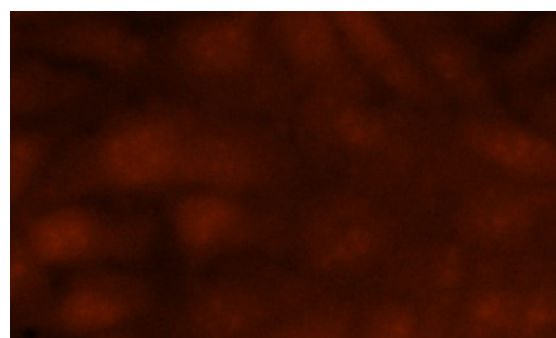

Rhod-2 AM

Figure 1. U2OS cells were seeded overnight at 40,000 cells per 100 µL per well in a 96-well black wall/clear bottom costar plate. The growth medium was removed, and the cells were incubated with 100 µL of 4 µM Rhod-2 AM or Rhod-4™ AM in HHBS at 37 °C, 5% CO<sub>2</sub> incubator for 1 hour. The cells were washed twice with 200 µL HHBS, then imaged with a fluorescence microscope (Olympus IX71) using TRITC channel.

## Directions for use

### Use of Rhod-4™ AM

#### Cell loading of Rhod-4™ AM esters

AM esters are the non-polar esters that readily cross live cell membranes, and rapidly hydrolyzed by cellular esterases inside live cells. AM esters are widely used for loading a variety of polar fluorescent probes into live cell non-invasively. However, cautions must be excised when AM esters are used since they are susceptible to hydrolysis,

[Info@fluoprobes.com](mailto:Info@fluoprobes.com)

[Technical-](mailto:Technical-support@fluoprobes.com)

[support@fluoprobes.com](mailto:support@fluoprobes.com)

[Order-online@fluoprobes.com](mailto:Order-online@fluoprobes.com)

Contact your local distributor

FluoProbes®, powered by



213 Avenue J.F. Kennedy - BP 1140  
93103 Montluçon Cedex - France  
Tél. 04 70 03 88 55 - Fax 04 70 03 82 60

P.2

FT-CQ6062

particularly in solution. They should be reconstituted just before use in high-quality, anhydrous dimethylsulfoxide (DMSO). DMSO stock solutions may be stored desiccated at -20°C and protected from light. Under these conditions, AM esters should be stable for several months.

Following is our recommended protocol for loading Rhod-4™ AM esters into live cells. This protocol only provides a guideline, should be modified according to your specific needs.

- Prepare a 2 to 5 mM stock solution of Rhod-4™ AM esters in high-quality, anhydrous DMSO. The nonionic detergent Pluronic® F-127 is sometimes used to increase the aqueous solubility of Rhod-4™ AM esters.  
*[Note: A 20% Pluronic F-127 solution can be used in replacing DMSO to prepare solutions of these calcium indicators. A variety of Pluronic F-127 solutions can be purchased from FluoProbes].*  
*[Caution: long-term storage of AM esters in the presence of Pluronic F-127 is not recommended].*
- On the day of the experiment, either dissolve Rhod-4™ solid in DMSO or thaw an aliquot of the indicator stock solution to room temperature. Prepare a working solution of 1 to 10 µM in Hanks and Hepes buffer (HHBS) or the buffer of your choice. For most of cell lines we recommend that 4-5 µM Rhod-4™ reagents be used. The exact concentration of indicator required for cell loading must be determined empirically. To avoid calcium buffering, toxicity and other artifacts of overloading, one should generally use the lowest probe concentration that yields sufficient signal.
- If your cells containing the organic anion-transporters, probenecid (1–2.5 mM) or sulfapyrazone (0.1–0.25mM) may be added to the cell medium to reduce leakage of the de-esterified indicators.
- Incubate cells with the Rhod-4™ AM esters for 30 minutes to one hour at room temperature or 37 °C.  
*[Note: Decreasing the loading temperature might reduce the indicator compartmentalization]*
- Wash cells twice in HHBS or buffer of your choice (containing an anion transporter inhibitor, such as 2.5 mM probenecid if applicable) to remove excess probes.
- Run the experiments at Ex/Em = 530/570 nm.

## Measuring Intracellular Calcium Responses

To determine either the free calcium concentration of a solution or the K<sub>d</sub> of a single-wavelength calcium indicator, the following equation is used:

$$[Ca]_{free} = K_d[F - F_{min}]/F_{max} - F]$$

where F is the fluorescence of the indicator at experimental calcium levels, F<sub>min</sub> is the fluorescence in the absence of calcium and F<sub>max</sub> is the fluorescence of the calcium-saturated probe.


The dissociation constant (K<sub>d</sub>) is a measure of the affinity of the probe for calcium. The Ca-binding and spectroscopic properties of fluorescent indicators vary quite significantly in cellular environments compared to calibration solutions. In situ response calibrations of intracellular indicators typically yield K<sub>d</sub> values significantly higher than in vitro determinations. In situ calibrations are performed by exposing loaded cells to controlled Ca<sup>2+</sup> buffers in the presence of ionophores such as A-23187, 4-bromo A-23187 and ionomycin. Alternatively, cell permeabilization agents such as digitonin or Triton® X-100 can be used to expose the indicator to the controlled Ca<sup>2+</sup> levels of the extracellular medium.

Trypan Red Plus™ can also be used to prevent Rhod-4 from leaking out of cells. It might inhibit the activities of drug-efflux pumps since it contains a probenecid-like moiety.

## Use of Rhod-4™ NW for HTS Applications

GPCR activation can be detected by direct measurement of the receptor mediated cAMP accumulation, or changes in intracellular Ca<sup>2+</sup> concentration. GPCR targets that couple via G<sub>q</sub> produce an increase in intracellular Ca<sup>2+</sup> that can be measured using a combination of Rhod-4™ reagents and a fluorescence microplate reader.

The fluorescence imaging plate readers has a cooled CCD camera imaging system which collects the signal from each well of a microplate (both 96 and 384-well) simultaneously. These plate readers can read at sub-second intervals, which enables the kinetics of the response to be captured, and has an integrated pipettor that may be programmed for successive liquid additions. Beside their robust applications for GPCR targets, our Rhod-4 Calcium Assay Kits can be also used for characterizing calcium ion channels and screening calcium ion channel-targeted compounds.

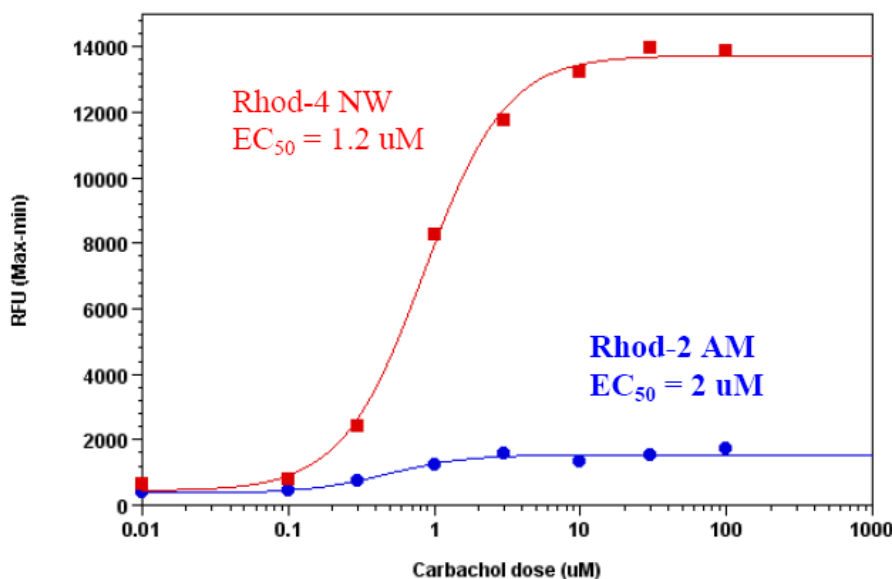

**Figure 2. Carbachol Dose Response in HEK-293 cells measured with Rhod-4 NW Assay kit and Rhod-2 AM under the same assay conditions.** HEK-293 cells were seeded overnight at 40,000 cells per 100  $\mu$ L per well in a 96-well black wall/clear bottom costar plate. The growth medium was removed, and the cells were incubated with 100  $\mu$ L of the Rhod-4 NW calcium assay kit, or Rhod-2 AM for 1 hour at room temperature. Carbachol (25  $\mu$ L/well) was added by NOVOstar (BMG LabTech) to achieve the final indicated concentrations. The EC<sub>50</sub> of Rhod-4 NW is about 1.2  $\mu$ M. The excitation and emission were 530nm and 570nm respectively.

Our Rhod-4 Calcium Assay Kits have the following advantages for HTS applications:

- Longer Wavelengths:* multiple excitations @ 488, 514, 532 & 546 nm; maximum emission @ ~555 nm
- No Wash Required and No Quencher Interference with Your Targets.*
- Robust Performance:* enable calcium assays that are impossible with Rhod-2 AM.
- Strongest Signal Intensity:* 4 times brighter than Rhod-2 AM.
- Larger Assay Window:* 10 times better than Rhod-2 AM

## Conclusions

Because of the importance of Ca<sup>2+</sup> in biology, numerous techniques/methods for analyzing the mechanisms of cellular and/or subcellular Ca<sup>2+</sup> activity have been established. Unfortunately, however, there is no one best technique/method with which one can measure Ca<sup>2+</sup>. Although each method for analyzing Ca<sup>2+</sup> activity has certain advantages over the others, each also suffers drawbacks. With the outstanding properties described above, we believe that Quest Rhod-4™ calcium detection reagents and Rhod-4NW Calcium Assay Kits provide new powerful tools for intracellular calcium analysis and monitoring in a variety of biological systems in coupling with the rapid advance in fluorescence instrumentation.

As might have been predicted, the interests of many researchers shifted from Ca<sup>2+</sup> analysis at the cellular level to that of the subcellular level. It has been found that Ca<sup>2+</sup> is not even distributed throughout the whole cell and that intracellular heterogeneity of Ca<sup>2+</sup> (such as Ca<sup>2+</sup> waves and Ca<sup>2+</sup> sparks) is observed in a variety of cells (e.g., oocyte, heart muscle cell, hepatocyte, and exocrine cell). With the advent of the confocal laser scanning microscope (CLSM) in the 1980s and advanced microplate readers dedicated for intracellular calcium detection) in 2000s, measurement of intracellular Ca<sup>2+</sup> has accelerated significantly. Confocal laser scanning microscopy, and more recently multiphoton microscopy, allows the precise spatial and temporal analysis of intracellular Ca<sup>2+</sup> activity at the subcellular level.

## Related products

- |   |   |
|---|---|
| - Trypan Red Plus, <a href="#">CL1090</a>   | - Rhod-4 NW, <a href="#">CQ6080</a>                       |
| - Fluo-3 AM, <a href="#">FP-78932C</a>      | - Fura-2 AM, <a href="#">FP-42776C</a>                    |
| - Fluo-8 NW for HTS, <a href="#">CJ2550</a> | - Pluronic F-127, 20% solution, <a href="#">FP-69806A</a> |
| - Indo-1 AM, <a href="#">FP-42775A</a>      | - Coelenterazine (native), <a href="#">UP972333</a>       |
| - Rhod-2 AM, <a href="#">FP-661583</a>      | - Coelenterazine 400a, <a href="#">UPBB8392</a>           |

[Info@fluoprobes.com](mailto:Info@fluoprobes.com)

[Technical-](#)

[support@fluoprobes.com](mailto:support@fluoprobes.com)

[Order-online@fluoprobes.com](#)

Contact your local distributor

FluoProbes®, powered by



213 Avenue J.F. Kennedy - BP 1140  
93103 Montluçon Cedex - France  
Tél. 04 70 03 88 55 - Fax 04 70 03 82 60

FT-CQ6062

- Coelenterazine H, [UPR3078B](#)
- A23187, [FP-28362A](#)
- 4-bromo A23187, [FP-37222A](#)
- Ionomycin, [FP-53989A](#)
- Probenecid, Cell culture tested, [FP-288652](#)
- Probenecid, water soluble, [FP-288653](#)
- Calcium calibration buffer kit, [FP-21527A](#)
- DMSO anhydrous, [JW7390](#)
- Black wall, clear bottom assay plate (96-well) [KT225](#), (384-well), [BA8170](#)

## Ordering information

Catalog size quantities and prices may be found at [www.interchim.com/](http://www.interchim.com/)

Please inquire for higher quantities (availability, shipment conditions).

For any information, please ask : FluoProbes® / Interchim; Hotline : +33(0)4 70 03 73 06

**Disclaimer :** Materials from ABD Bioquest are sold **for research use only**, and are not intended for food, drug, household, or cosmetic use.

**Warning:** This kit is only sold for the end users. It is covered by a pending patent. Neither resale nor transfer to a third party is allowed without written permission from ABD Bioquest. Chemical analysis of kit components is strictly prohibited.

[Info@fluoprobes.com](mailto:Info@fluoprobes.com)

[Technical-](#)

[support@fluoprobes.com](mailto:support@fluoprobes.com)

[Order-online@fluoprobes.com](http://Order-online@fluoprobes.com)

Contact your local distributor

FluoProbes®, powered by



213 Avenue J.F. Kennedy - BP 1140  
03103 Montluçon Cedex - France  
Tél. 04 70 03 88 55 - Fax 04 70 03 82 60

P.5