




Instructions are shown using  
item no. 45237 as an example

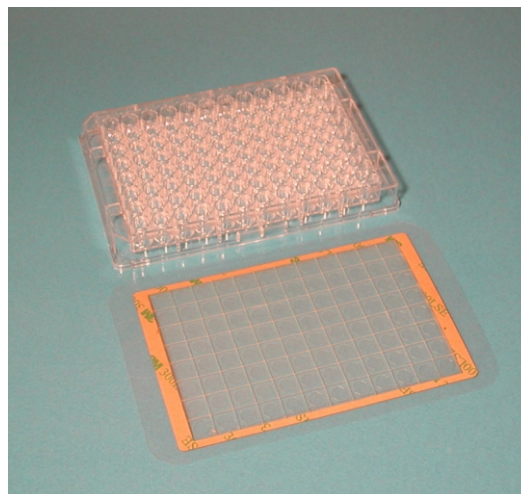


1. Plate and Cover  
Cover: adhesive liner facing up.

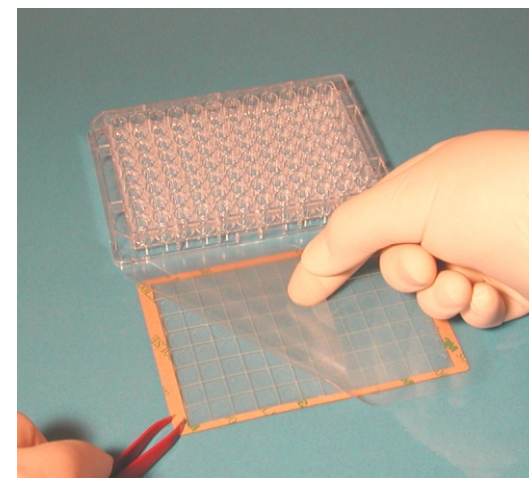

2. To expose the adhesive; peel off  
the clear release liner.

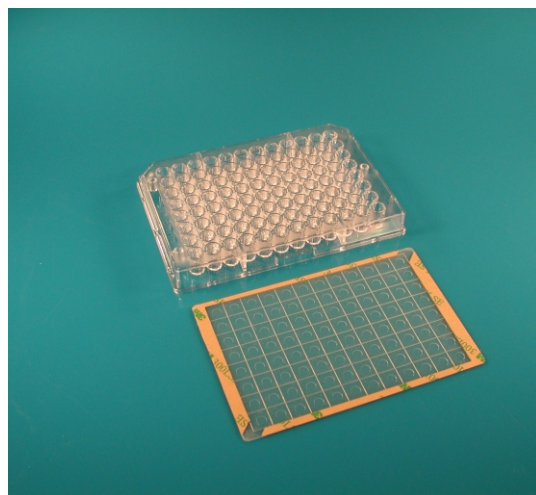

3. Release liner completely removed

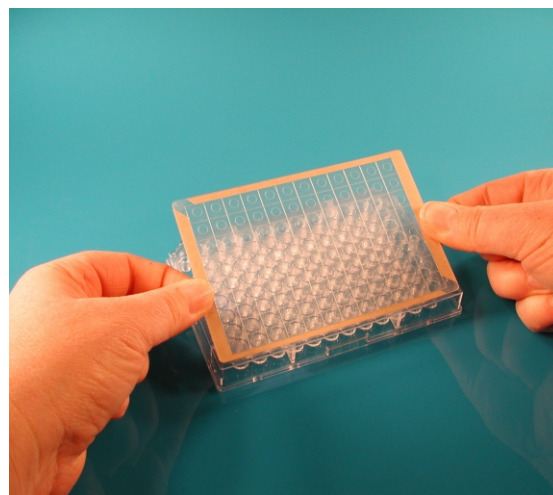

4. Handle cover by the protected edges,  
aligning to the plate over the wells.




5. Press firmly to seal cover to the  
wells.


Grace Bio-Labs, Inc.

Tel: 541-318-1208 / Fax: 541-318-0242 / E-mail: [custservice@gracebio.com](mailto:custservice@gracebio.com) / [www.gracebio.com](http://www.gracebio.com)




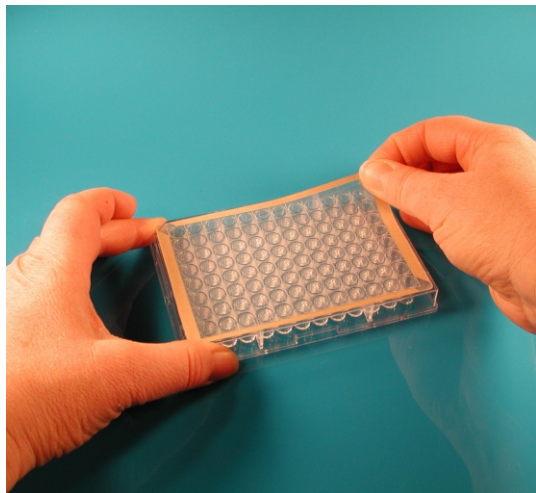

6. Remove backing liner and frame by peeling slowly.

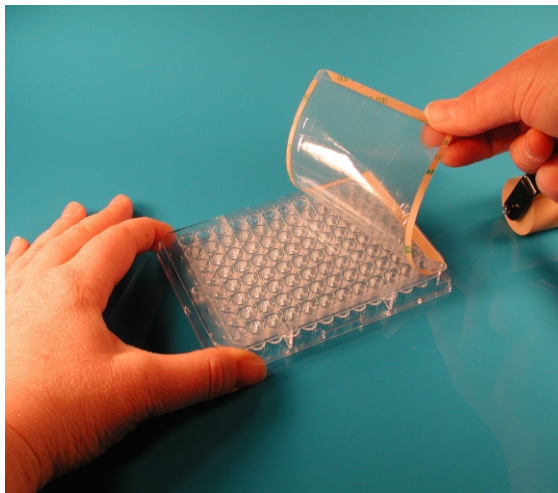

7.

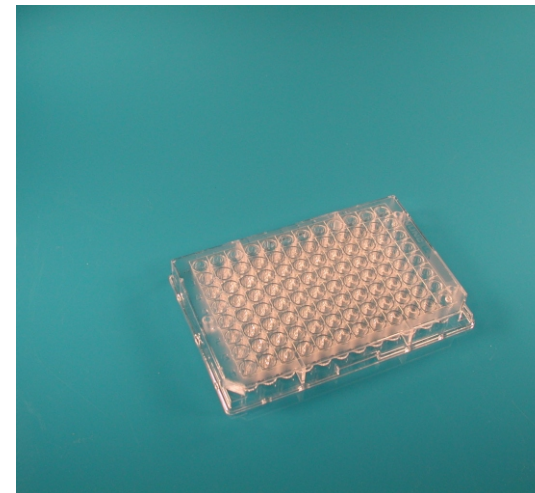


8. Frame & backing liner completely removed.



9. Remove an individual well cover by inserting forceps in the corner, grasp firmly and peel off.



10.

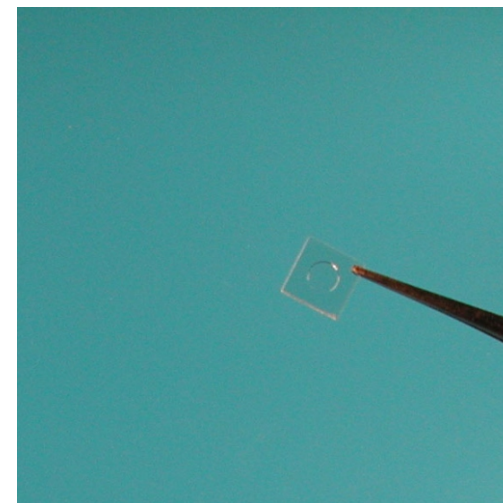

11.

Grace Bio-Labs, Inc.

Tel: 541-318-1208 / Fax: 541-318-0242 / E-mail: [custservice@gracebio.com](mailto:custservice@gracebio.com) / [www.gracebio.com](http://www.gracebio.com)

