

## CultureWell™ Gaskets







### Product Information

Removable Chambered Coverglass, non-sterile

Cat. number, Qty	Silicone Well Depth (mm)	Volume Per Well (µl)	
FP-JZ7603, 10 U	1.0	250 - 400	
FP-CK6460, 10 U	1.0	300 - 500	
FP-CK6470, 10 U	1.0	50 - 100	
FP-CK6480, 10 U	1.0	15 - 30	
FP-CK6490, 10 U	1.0	3 - 10	

## Directions for use

1. Cell culture may be performed using conventional methods.
2. Gently decant media.
3. Separate the two black gasket tabs and insert the slip separator string between gasket tabs as shown.
4. Pull slip separator along the length of the assembly to separate silicone coverglass gasket and chambered well gasket.
5. Gently remove coverglass with attached gasket.
6. Small volume incubation may be performed using the wells formed by the silicone gasket.
7. Aqueous mounting of specimens may be performed with wells in place, by filling wells with media and pressing a coverglass or microscope slide to the gasket surface.
8. To remove the silicone gasket from the coverglass, peel slowly at a sharp angle to the glass.
9. For specimen mounting, match the less rounded corner of the coverglass to location feature (H) on the blue frosted edge of the microscope slide (provided). Refer to orientation key below.

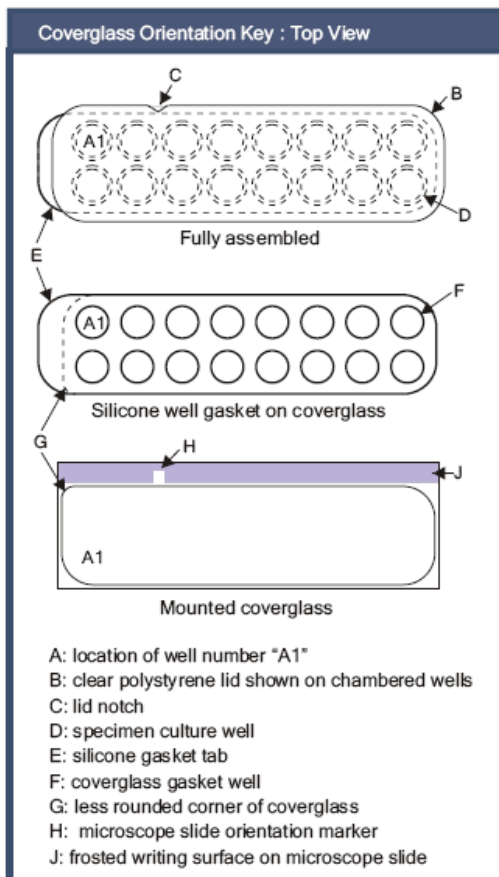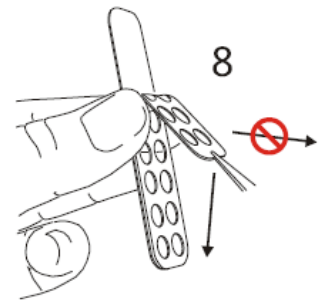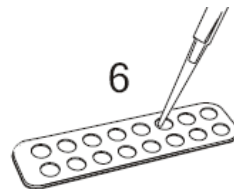
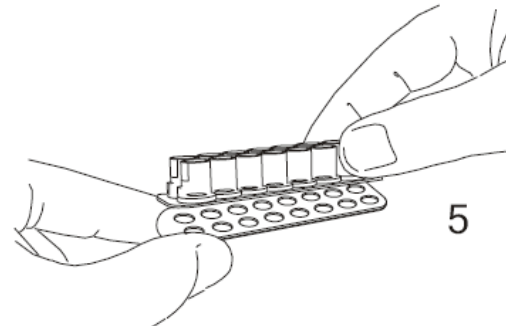
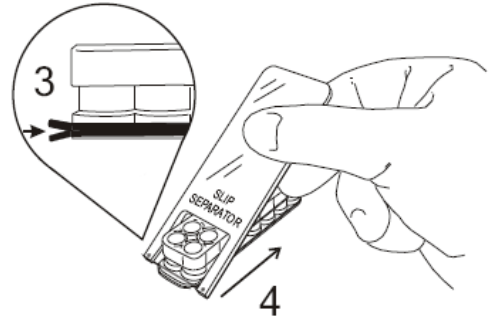
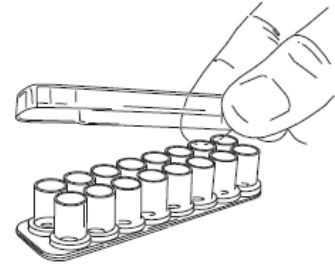


FT-JZ7607

CultureWell™  
Removable Chambered Coverglass  
Data Sheet

EXPERIMENT NO.:

DATE:

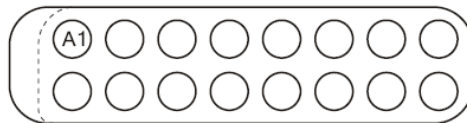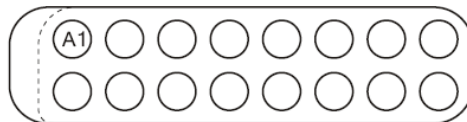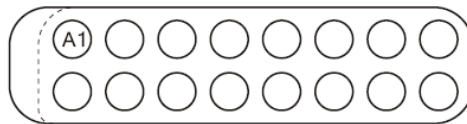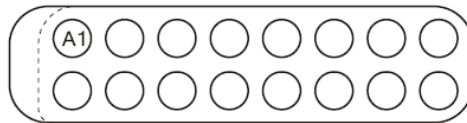

## Ordering information

Catalog size quantities and prices may be found at <http://www.interchim.com>  
Please inquire for higher quantities (availability, shipment conditions).

For any information, please ask : FluoProbes® / Interchim; Hotline : +33(0)4 70 03 73 06

**Disclaimer :** Materials from FluoProbes® are sold **for research use only**, and are not intended for food, drug, household, or cosmetic use.  
FluoProbes® is not liable for any damage resulting from handling or contact with this product.

[Info@fluoprobes.com](mailto:Info@fluoprobes.com)  
[Technical-support@fluoprobes.com](mailto:Technical-support@fluoprobes.com)  
[Order-online@fluoprobes.com](mailto:Order-online@fluoprobes.com)

Contact your local distributor

FluoProbes®, powered by



213 Avenue J.F. Kennedy - BP 1140  
03103 Montluçon Cedex - France  
Tél. 04 70 03 88 55 - Fax 04 70 03 82 60